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**W16**, a water bath for temperatures up to 110°C. The container, made of special steel, is compactly built into a container which is also constructed from special steel. The temperature is controlled by a Thermostat which is exact up to +- 1,5 K and more.

A dry-go protection prevents overheating when there is a water shortage. Added protection is also available with the temperature fuse. A deep lying Ring set (Order No. RI 17/E) allows the condensed water to fall back into the bath.

**PZ 26-4**, which a capacity of 5 litres, can be used as fluid, air and sand bath up to 400°C. As a bath-filling, the combinations of oils, watersoldering tins, sand or other bath-fillings can be used without detrimentally affecting the bath, wh@ is made out of special steel - W.No. 1.4301.

**PZ 26-7**, designed for temperatures of up to 700°C offers the same possibilities but not, however, as a fluid bath. This apparatus can only be used as a dry air or sand bath. Both baths are also well-suited for use as rotary evaporators.

PZ26-4 and PZ26-7 are both heated with a tubular radiator.

**MB 28/P** is a fluid bath with an especially large immersion depth of 280 mm, with a bath diameter of 220 mm. The special steel bath container, material No. 1.4301 has no bore holes, a large radius at the bottom and is equipped with a 6-step switchable underfloor heating device switchable performances: 135-220-300-850-1150-1500 Watt. The instrument is equipped with a built-in-thermo-monitor.

PZ26.. and MB28/P are microprocessor controlled by a PID Regulator with a low-mass sensor. The actual value temperature is allways on view. To see the temperature setpoint a button has to be pressed. The setpoint temperature is adjustable with an up/down-button.

The housing is constructed from special steel and the base is lacquered. All types of bath are constructed and build for continuous operation, even at maximum temperatures.

Their housings and all component parts will heat up only slightly.

W16, P26-4 u. PZ26-7 in equipped with a screw-on fixture for a support rod of 12 mm dia. MB 28/P for 2 support rod.

Technische Daten Best.-Nr. = Typ		<b>W 16</b>	<b>PZ 26-4</b>	<b>PZ 26-7</b>	<b>MB 28/P</b>	Technical Data Order No. = types
Temperaturvorwahl	°C	30...110	20...400	20...700	20...240	Range of temperature
Anzeige		ohne/without	digital	digital	digital	Display
Leistung	W	1000	1200	1200	1500	Performance
Spannung	V	220...240 V, 50-60 Hz				Voltage
Außenabmessung	mm	220 x 220	310 x 310			Overall dim.
Höhe	mm	160	280	280	435	Height
Gewicht	kg	3,0	6,8	6,8	10,5	Weight

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### Further Fabrications

High temperature TITANIUM hot plate up to 600°C - Precision hot plates - Stand hot plates with supporting rod - rapid incinerator - Multi-purpose, extremely deep, heating bath - sand bath - Series heating units - Multipurpose heating equipment - Water- and Oil baths - Development - Special orders.

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The right to make modification serving further technical development, is reserved