AGITADOR ROTATIVO «MOVIL ROD»

TUBE SHAKER «MOVIL ROD»

7001723

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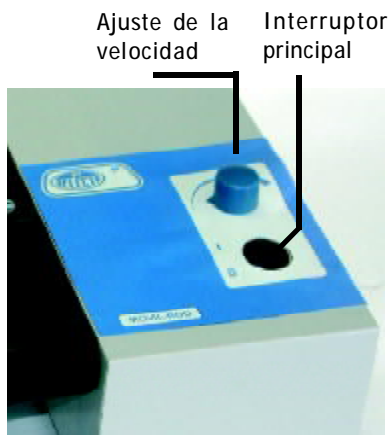
2 Installation

Unit description:

Shaker with 7 rollers with double motion action: rotation and swing, this provide a suitable homogenize effect.

Suitable for use in liquid suspension, biology and haematology

Application examples:

- Suitable for mixing blood and anticoagulant.
- Mixing of substances as dry powder or frozen.
- Liquid mixing.

Location:

- Place on flat and levelled surface. Suitable for unit weight.
- Let free 20cm, at least, around the unit.
- Near to power socket (1m).

Connection to power supply:

- Check that power socket have suitable voltage, according unit plate.
- Power socket have earth connector. Unit must be grounded.
- Unit could be dangerous is not properly grounded.

3 Technical specification

Rolling and vibrating speed:	From 10..80 rpm.
Vibrating range:	22mm±1mm
Power consumption:	12W
Fuse:	2A (Ø5 x 20mm)
Voltage:	230V AC ±10% (50/60Hz)
Weight:	3.6Kg
Dimensions:	394 x 266 x 98 mm

4 Operation

- Place the tubes to stirr.
- Press the main switch
- Adjust the speed.

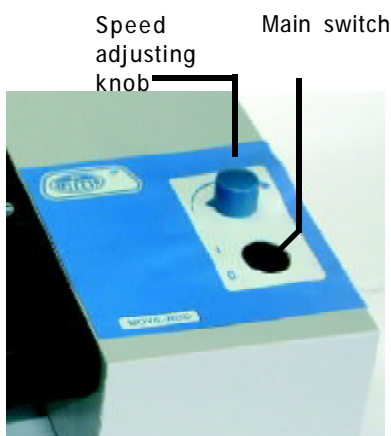
Near once a year, lubricate ,slightly, the bearings.

5 Warranty

The warranty covers the defective components or faulty assembling of the unit for 1 year since the purchasing date. Keep purchasing bill as the only valid document for the purchasing date.

The warranty does not covers the damages caused for a unit malfunction.

Return goods are not accepted without previous J.P. SELECTA,s.a. approval.

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Electrical equipment for measurement, control and laboratory use. EMC Requirements.

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Safety requirements for electrical equipment for measurement, control and laboratory use. Part 1: General requirements.

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Particular requirements for laboratory equipment for mechanical mixing and stirring.

Abrera, Enero 2007

Ramón Ramón
Dtor. Técnico.

David Pecanins
Rble. Calidad