## CENTRÍFUGAS - CENTRIFUGES

**CENTRONIC BLT**

**7001768**

**CENTROFRIGER BLT**

**7001769**

**CENTROMIX BLT**

**7002556**

**MIXTASEL BLT**

**7002576**

**MEDITRONIC BLT**

**7002774**

# Indice

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## 1 Safety

The equipment includes appropriate safety measures. Risky situations that must be respected are indicated along this manual.

### 1.1 Safety icons

The safety icons identify the risky situations and the safety measures that must be taken.

Risk of danger  
Electrical risk

Respect the indicated instructions to perform the operation.

### Electrical risks

Risk of electrical accident when accessing the areas indicated by this sign or when performing the operations indicated by this icon in this manual.

Respect the indicated instructions to perform the operation.

### Important information

- Important information for good results or for an optimum equipment operation.
- Important information to extend life of the equipment or to prevent degradation of any of the components.

### 1.2 Personnel qualification

This equipment can only be used by personnel who have been properly qualified.

This equipment can only be used by personnel who have read and understood these instructions or who have been properly qualified in the operation of this equipment.

## 2 General information

Handle the package with care. Unpack and check that the contents coincides with the "Packing List". If any part is damaged or missing, please advise the distributor immediately.

Do not install or use the equipment without previously reading this instructions manual. It is an inseparable part of the equipment and should be available to all users of the machine. Any questions can be clarified by contacting J.P. SELECTA's technical service.

Any modification, deletion or lack of maintenance of the device could cause damages. The manufacturer is not liable for the damages that may arise.

Do not use the equipment with fluids which can give off vapours or form explosive or flammable mixtures.

### Maintenance

**Manipulation of electronic circuits by unauthorized persons may cause damages difficult to repair. Check with J.P. SELECTA's technical service.**

To clean the accessible parts, use cleaner or a similar product. Do not use aggressive solvents.

### Guarantee

This product is guaranteed for one year. The warranty does not cover damages caused by an improper use or causes beyond J. P. SELECTA, s.a.u.

Any manipulation of the equipment by an unauthorized personnel cancels the warranty benefits.

### 3 Technical specifications

**Voltage supply 115/230V 50/60 Hz as indicated in the machine's features plate.**

Code:	7001768	7001769	7002556	7002576
Dimensions (mm): Height x Width x Depth	340x410x540	340x410x710	230x300x350	340x390x440
Power (W):	440	960	180	210
Max speed (rpm):	13500	15300	4400	4200
Max RCF: (Depending on the rotor)	19048	23700	2791	3151
Weight (kg):	22	39	12	22
				25

### 4 Packing List


#### 4.1 Standard package content

The standard equipment consists of the following components:

- Centrifuge CENTRONIC BLT: 7001768
- Centrifuge CENTROFRIGER BLT: 7001769
- Centrifuge CENTROMIX BLT: 7002556
- Centrifuge MIXTASEL BLT: 7002576
- Centrifuge MEDITRONIC BLT: 7002774
- Network connection cable: 7140
- Rotor nut wrench\*: 5377
- (\*only for models: CENTRONIC BLT  
CENTROFRIGER BLT  
MIXTASEL BLT)
- Instructions manual 80315

#### Notice to customers: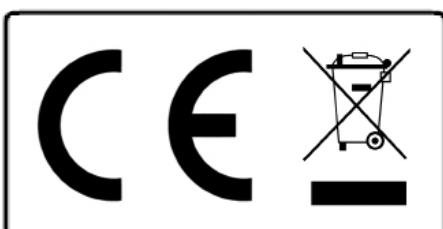

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

## 5 Introduction

The centrifuge has an induction motor maintenance free and a maximum noise level of 60 dBA. It has an internal security camera and a bucket, stainless steel cover.

The electronic circuit is controlled by a microprocessor that controls the operating parameters: speed, acceleration, braking, RCF, temperature and operating time. It can store up to 10 operating programs.

FREE SPACE AROUND CENTRIFUGE

### Safety zone:

Leave a free space of 30 cm. around the centrifuge.

With the centrifuge running:

- Do not place hazardous substances in the security area.
- The staff should be outside the safety area.

## 6 Installation

### 6.1 Location

- Place the equipment near a power socket suitable for consumption and over a flat level surface and strong enough to support the equipment weight and possible vibrations.
- The surface must allow the adherence of the centrifuge legs.
- Leave a free space of 30 cm. around it (See danger indication).
- This equipment can cause vibrations that could affect other nearby equipments.
- Do not install the centrifuge near heat sources or in places where it is exposed to direct sunlight, as this may affect temperature regulation.
- Place the nipple provided in outlet (16) and connect to a hose to drain condensates.

**DO NOT INSTALL THE EQUIPMENT IN EXPLOSIVE ATMOSPHERES OR PLACES WHERE SUBSTANCES LIABLE TO FORM EXPLOSIVE OR FLAMMABLE MIXTURES ARE USED OR STORED.**

### 6.2 First lid opening and emergency opening

Upon receiving the centrifuge, the network connection cable is inside the tank and therefore the only way to open the equipment is following these instructions:

- Remove the equipment side plastic caps.
- Gently push from both caps and lid will open.

In MIXTASEL BLT models, remove the plastic cap located at the bottom of the equipment. Pull gently on the cap and the lid will pop open.

**USE THIS OPENING SYSTEM FOR EMERGENCY OPENING ONLY.**

**NEVER use the device emergency opening when the engine is rotating.**

### 6.3 Connection to the mains

Make sure that the equipment is connected to the mains and that the voltage coincides with the indicated on the features plate.

Do not use an equipment without being earthed.

Only use the supplied cable for network connection.

### 6.4 Rotor positioning

Before starting the centrifuge, ensure that you have placed the rotor in a proper way, otherwise it could cause damages on the equipment.

Make sure to fit the pin of the motor shaft in its housing.

Firmly tighten the nut of the rotor with the provided key (*only for models Centronic BLT, Centrofriger BLT and Mixtasel BLT*).

Place the adapters and tubes evenly.

## 7 Operation

All the functions of this equipment are set by the touch screen panel. Softly press with the finger in the centre of each icon, and the action will be performed when releasing the finger from the screen.

When starting the equipment, the starting screen is displayed.

- **Settings:**


- Time and date: menu to set date and time.

- **Programming:**


- Enables programming the centrifuge profile, by defining speed, time and acceleration and deceleration ramps.

- **Task:**

- Enables selection and implement a centrifuge profile.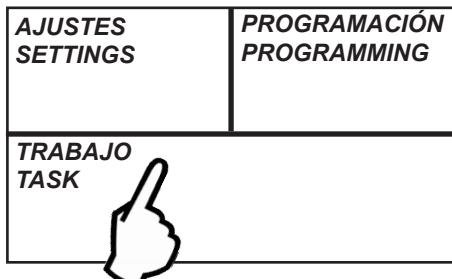


## Screen diagrams:

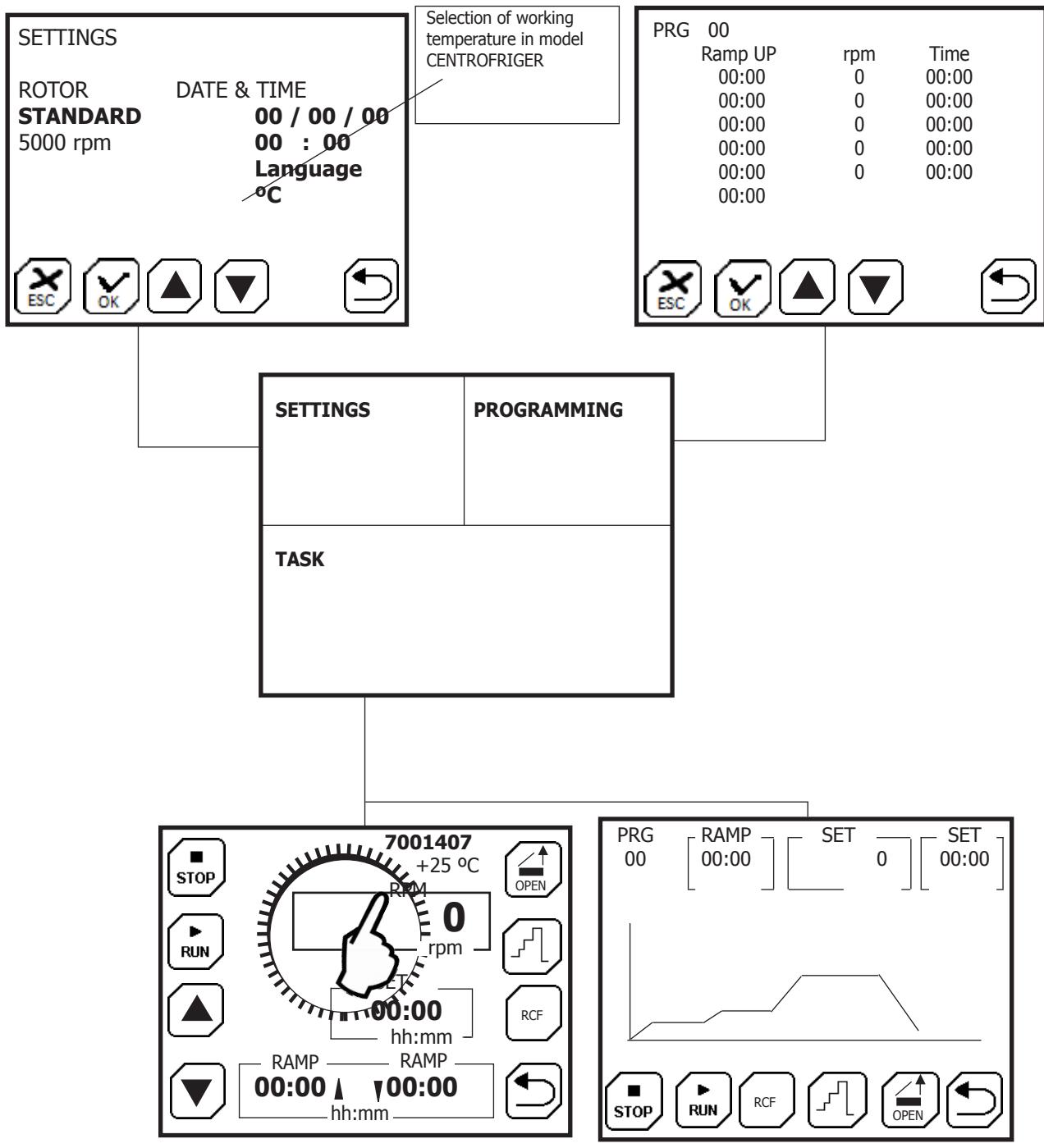To change mode, press
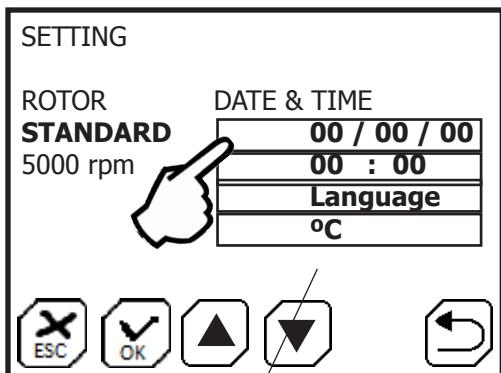
### 7.1 Setting screen

- Update date, time.  
Press on the Date and Time panel.

Use the keys:


and accept by pressing:

In a similar way:

- Select the Language.
- Select the Temperature units.

Selection of working temperature in model CENTROFRIGER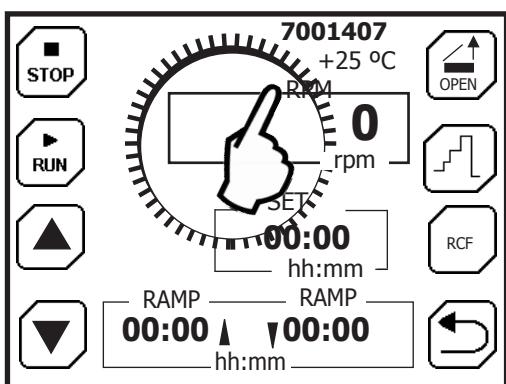

### 7.2 Direct programming and operation


- Press on the RPM display.
- Press on to set the RPM.
- Press on SET to establish minutes/hours of operation.
- If necessary, set the acceleration and deceleration times of «RAMP» in
- RAMP: Acceleration time up to rpm set point.
  - RAMP: Deceleration time.

(Setting ramp minutes) A time of 00.00 indicates the fastest.

Press to start centrifugation

The same screen indicates the centrifugation process evolution.


Once centrifugation is finished, open the lid with: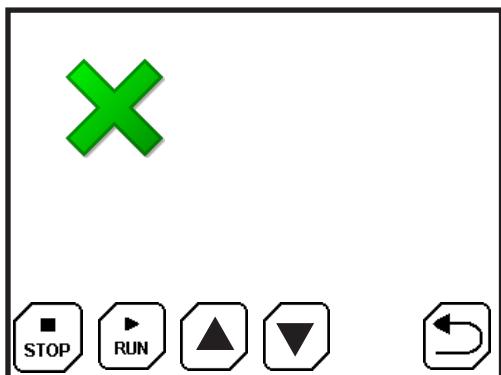

### 7.3 Error message

Any incidence in operation causes an error message.

Fig 7.3 Error screen

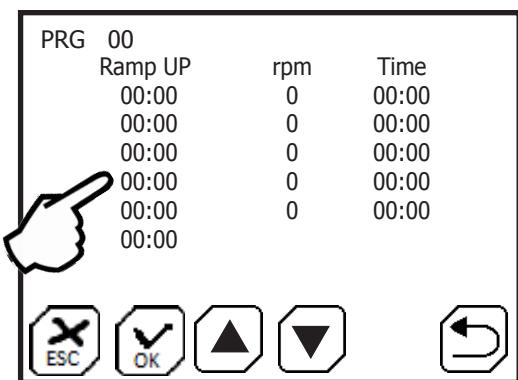## 7.4 Profile programming

To program a profile, press:

A speed profile is formed by two segments.

One segment establishes:

1. **RAMP time** to reach speed.
2. Centrifugation **speed** (rpm).
3. Centrifugation **time** (hh:mm).

If RAMP time is zero or very small, the centrifuge will be conducted in the quickest way.



A segment with time 0 is not executed.

The last ramp time is braking to stop the rotor.

### Centrifugation profile example

PRG 01

Ramp UP	rpm	Time
00:10	1000	00:25
00:15	2000	01:15
00:35		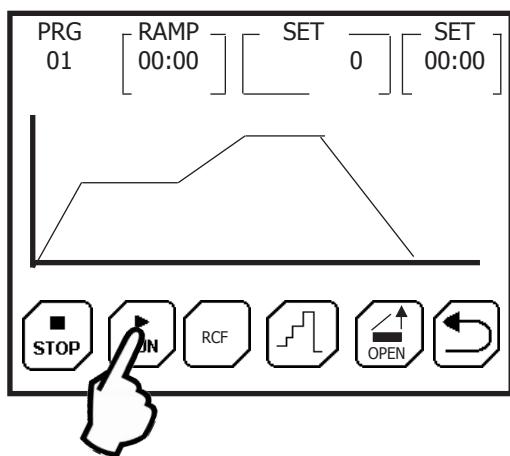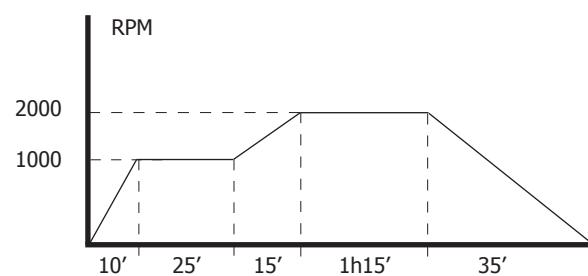

The process monitoring screen indicates at any time the ramp time, speed (rpm) and centrifugation time.