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**ELECTRONIC-DIGITAL CONTROL CENTRIFUGES
WITH INDUCTION MOTOR**

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Index

1 Safety	10
1.1 Safety Icons	10
1.2 Risks that the operator can be subjected to	10
1.3 Qualified users	10
2 General Information	11
3 Technical Specifications	11
4 Contents List	11
4.1 Standard contents	11
5 Introduction	12
6 Equipment Description	12
6.1 Equipment	12
6.2 Features	12
6.3 Control panel	12
6.4 Rear part	12
7 Installation	13
8 Operation	13
9 Maintenance	14
10 Guarantee	15
11 Electric Diagram	16

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1 Safety

The equipment comes complete with the safety features.

This manual indicates areas of possible risk.

1.1 Safety Icons

Identifies risk situations and safety measures that should be adhered to. The icons make reference within the paragraph, marked as a lined grey.

Risk danger

Risk danger

Beware, follow instructions as indicated.

Electrical risk

Risk of electrical shock if accessing zones are shown with this icon.

Beware, follow instructions as indicated.

Important information

- Important information in obtaining the best results or optimum performance of the equipment.
- Important information to extend the life of the equipment and maintain its optimum performance.

1.2 Risks that the operator can be subjected to

- Blows risk.
- Electrical risk.

1.3 Qualified users

This equipment should only be used by personnel that are qualified users.

This equipment should only be used by personnel that have understood this manual or have previous adequate instruction of the use of centrifuges.

2 General Information

- 2.1 Handle the equipment with care. Unpack and check that all items coincide with the delivery note. If you see any discrepancy or damage of any kind, notify your distributor as soon as possible.
- 2.2 Read this manual before using the equipment.
- 2.3 These instructions are an important part of the equipment and should be made available to all users.
- 2.4 If in any doubt or clarity on the use of this instrument, contact your local distributor or the service department of J.P. Selecta, s.a.u.
- 2.5 **ATTENTION, EQUIPMENT NOT CLEAN AND DISINFECTED SHALL NOT BE REPAIRED.**
- 2.6 Any unauthorised modification, lack of maintenance or deliberate damage to the machine, transgresses the directive 89/655/CEE where the manufacturer will not be responsible for damages that could occur due to undue care and attention.
- 2.7 Do not use the equipment for fluids that can produce vapours that can form inflammable or explosive mixtures.

3 Technical Specifications

Voltage supply 115-230V 50/60Hz according to the characteristics plate indications.

- Code: 7002356
- Dimensiones (cm):
 - High: 20
 - Width: 28
 - Depth: 34
- Power (W): 170
- Speed max. (rpm): 5000
- RCF max.: 2515
- Weight (kg): 6

4 Contents List

4.1 Standard contents

The standard equipment consists of the following items:

- Centrifuge: 7002356
- Angle head for 8 tubes 15ml 30098
- 8 Aluminium buckets of 15ml 7001500
- 8 Glass tubes of 15ml 7000772
- Power supply cord 7140
- Instruction manual 80224

5 Introduction

The centrifuge Centro-8-BL is a tabletop centrifuge with a free maintenance induction motor and electronic digital control of speed and working time, controlled acceleration and deceleration, temporized mode from 1 minute to 60 minutes and safety electro-mechanic lock.

6 Equipment Description


6.1 Equipment

The Centro-8- BL centrifuges have the following features:

- Electronic digital control circuit board
- Induction motor (brushless)
- Angle rotor 8 tubes of 15 ml
- Electromagnetic safety lock

6.2 Features


- Regulable speed until 5000 rpm
- Max RCF 2515 g
- Speed reading on the display (resolution 10 rpm)
- Working time reading on display (resolution 1 min)
- Instant start/stop key to the set speed
- Electronic lid opening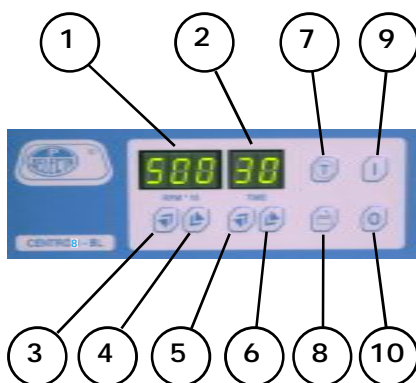

6.3 Control panel

1. Speed display (rpm)
2. Time display (min)
3. Increase speed value key
4. Decrease speed value key
5. Increase time value key
6. Decrease time value key
7. Instant start/stop key
8. Start key
9. Stop key

6.4 Rear part

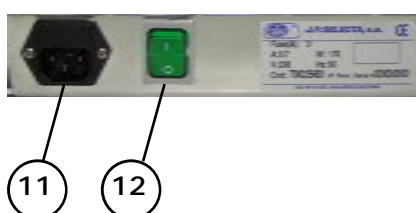10. Power supply socket
11. Main switch

7 Installation

7.1 PLACEMENT

Place the piece of the equipment near of a socket, keeping in mind its amperage, on a leveled surface enough resistant to support the weight indicated on the table of the section 3 (Technical features), leaving and space of 10cm by the rear part and both laterals. If it is a table, make sure that its legs are able to absorb the centrifuge vibrations.

DO NOT INSTALL THE CENTRIFUGE IN AREAS WHERE INFLAMMABLE LIQUIDS ARE STORED OR SPECIAL PROTECTION AREAS.

7.2 POWER SUPPLY CONEXION

Make sure that the voltage given by the socket is the indicated on the feature label attached on the rear part.

Do not use the equipment unless it is correctly earthed.

If you replace the plug keep in mind the following:


Blue wire:	Neutral
Brown wire:	Phase
Yellow/green wire:	Ground

8 Operation

STARTING UP

- Connect the equipment to the power supply
- Turn on the main switch (12). The displays (1) and (2) will light.
- Press key (8) to open the lid.
- Insert the tube in the rotor in a balanced compensate way.
If only use 2 tubes, insert them in oposite sides of the rotor (not contiguous)

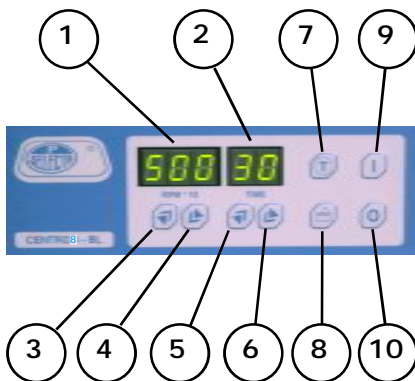**Never place one or three tubes.
Always two or four tubes.**

- Close the lid pressing on it until you heard the sound of the lock.
- Pressing the keys (3) and (4) increases or decreases the speed set which is shown on the display (1).
- Pressing the keys (5) and (6) increases or decreases the time set which is shown on the display (2).

WORKING FOR A NOT DEFINED TIME

When the display (2) shows «00», by pressing the key (6) the display shows: «--». If the centrifuge starts with this set time it only will stop manually.

- Start pressing the key (7). The centrifuge will stop once finished the time set.
During the operation, the display (1) shows the speed read by the tachometer.
The operation can be stopped by pressing the key (10).

INSTANT START/STOP

- Do the previous operation as explained on the section «Starting up». Once the speed and time of the operation are set and they are different of 0, by pressing the key (7) the rotor starts to turn at the speed set. By stop pressing, the centrifuge will stop.

EMERGENCY OPENING

In case of a cut of the power supply and it would be necessary to open the door, do as follows:

- Remove the plastic stopper on the right lateral of the equipment.
- Pull carefully from this stopper and the lid will open.

Do not use this emergency opening device when the machine is operating.

ERRORS LIST:

- Error 1: Eprom failure
- Error 3: Tachometer speed failure
- Error 4: Overcurrent
- Error 5: Low voltage
- Error 6: Motor feedback voltage too high

9 Maintenance

Before removing the cover to handle inside, unplug the equipment from the socket.

The handle of the electronic circuit boards by unauthorized personnel may cause damages. Make sure to give the equipment to an authorised technical assistance.

9.1 CLEANING

- Regularly clean the centrifuge chamber and, if necessary, clean with soap or a mild detergent and a damp cloth. This serves to hygiene and at the same time prevents corrosion by accumulations of dirt.
- Use commercially-available disinfectants such as, for example, Sagrotan, Buraton, or Terralin (available at chemist's shops or drugstores).
- The centrifuge and the accessories consist of various materials. A possible incompatibility must be considered.
- Before using cleaning or decontamination agents that were not recommended by us, contact the manufacturer to ensure that such a procedure will not damage the centrifuge.
- For autoclaving, consider the continuous heat resistance of the individual materials (see "Autoclaving" below).

Autoclaving

The service life of the accessories essentially depends on the frequency of autoclaving and use.

- Replace the accessories immediately when the parts show changes in colour or structure or in the occurrence of leaks etc.
- During autoclaving, the caps of the tubes must not be screwed on in order to avoid the deformation of the tubes.

Accessories	Max Temp.	Min Time	Max Time	Max cycles
Aluminium buckets	134-138	3	5	-
Aluminium rotors	134-138	3	5	-
Glass tubes	134-138	3	40	-
Polyallomer / polycarbonate rectangular carriers	115-118	30	40	-
Polyallomer / polycarbonate round carriers	115-118	30	40	-
Polyamide buckets	115-118	30	40	10
Polycarbonate / polyallomer lids for angle rotors	115-118	30	40	20
Polycarbonate tubes	115-118	30	40	20
Polyphenylsulfone (PPSU) caps for buckets	134-138	3	5	100
Polypropylene balance adapter for blood-bag systems	115-118	30	40	n.s
Polypropylene copolymer tubes	115-121	30	40	20
Polypropylene rectangular carriers	115-118	30	40	-
Polypropylene rotors	115-118	30	40	20
Polypropylene round carriers	115-118	30	40	-
Polysulfone caps for buckets	134-138	3	5	100
Polysulfone lids for angle rotors	134-138	3	5	100
Rubber adapter	115-118	30	40	-
Stainless-steel balance weight for blood-bag systems	121	30	30	n.s
Teflon tubes	134-138	3	5	100

9.2 SPARE PARTS

The more commons spare parts are:


Spare description	7002356
Electromagnetic lock	10040
Main switch	20017
Motor	24226
Tachometer sensor circuit board	29230
Control panel circuit board	29364
Power circuit board	29366
Mains filter	45011
Transformer 115-115/18V 1.2A	45048

10 Guarantee

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, s.a.u.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, s.a.u. automatically cancels the guarantee.

11 Electric Diagram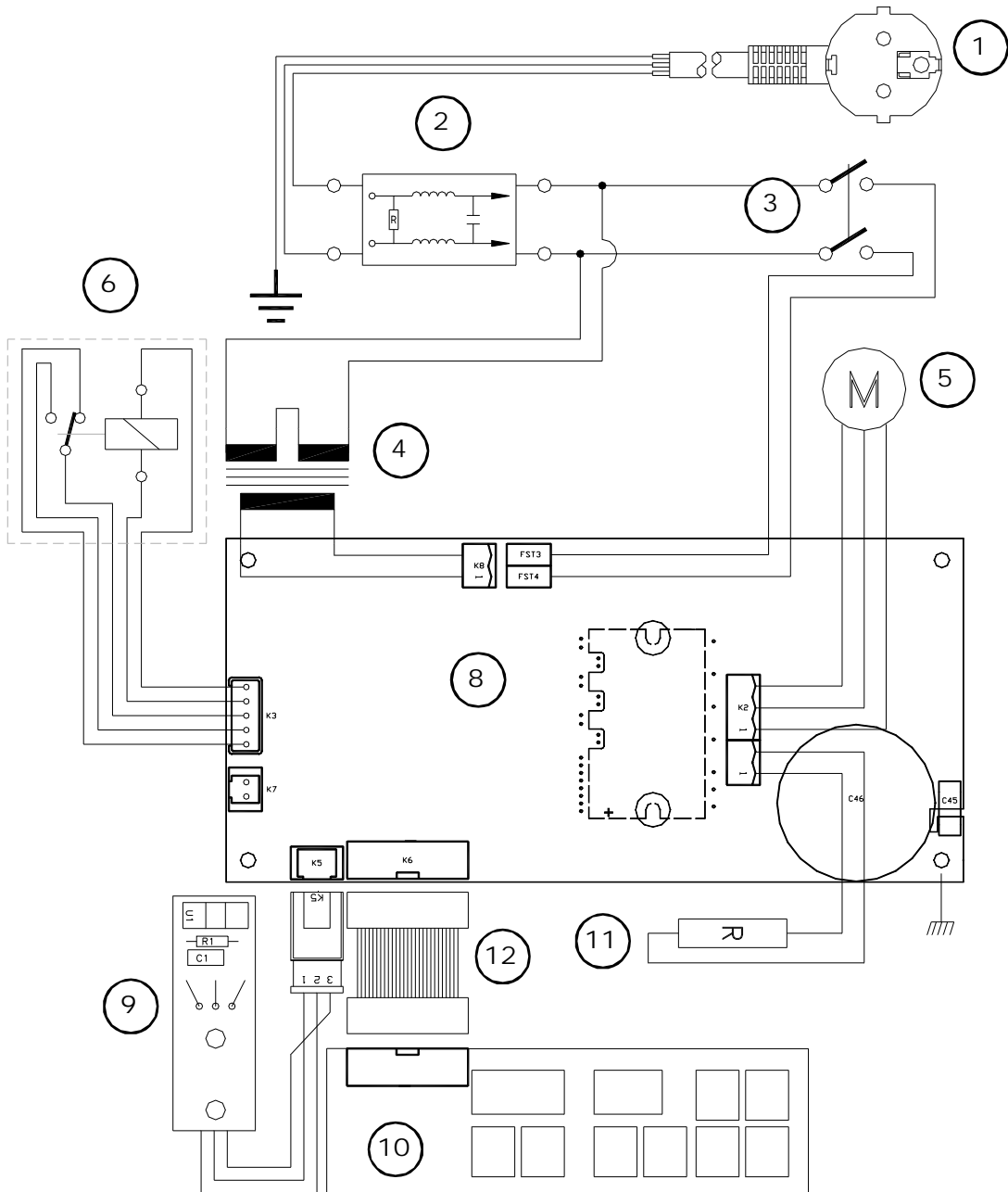

Pos	Description	Code
1	Power supply cord	7140
2	Mains filter	45011
3	Main switch	29017
4	Transformer 115-115/18V 1.2 A	45048
5	Motor	24226
6	Electromagnetic lock	10040

Pos	Description	Code
8	Power circuit board	29366
9	Tachometer sensor circuit board	29230
10	Control panel circuit board	29364
11	Break resistance	
12	Board connecting wire	