DRYTERM

2000787

**ESTERILIZADOR POUPIEL CALOR SECO
POUPINEL STERILIZER**

INFORMACIÓN GENERAL

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY EQUIPMENT TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the equipment with liquids which can give off vapours capable of making explosive mixtures.*

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The standard equipment consists of the following components:

Pieza	Piece	Código / Code	Cant. / Qty.
Esterilizador Dryterm	<i>Dryterm</i>	2000787	1
Bandejas	<i>Shelves</i>		2
Manual de instrucciones	<i>Instructions Manual</i>	80026	1

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*Poupinel sterilizers **DRYTERM** are designed for quick sterilization of diverse instruments (surgical, odontological, etc...) for adjustable temperatures from 60°C up to 250°C with a temperature fluctuation of $\pm 6^\circ\text{C}$.*

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Heating by shielded elements in the base which provide a quick temperature rise.

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Voltage supply 115-230 V 50/60 Hz as indicated on the characteristics plate.

Código / Code	2000201	
Capacidad / Capacity (L)	19	
Temperatura máxima / Maximum temperature	250°C	
Homogeneidad* / Homogeneity*	$\pm 4^\circ\text{C}$	
Estabilidad* / Stability*	$\pm 10^\circ\text{C}$	
Medidas interiores (cm) / Internal dimensions (cm)	Alto / Height	25
	Ancho / Length	32
	Fondo / Width	23
Medidas exteriores (cm) / Overall dimensions (cm)	Alto / Height	37
	Ancho / Length	54
	Fondo / Width	34
Consumo / Consumption (W)	770	
Peso / Weight Kg.	19	

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* *Measured according to the measurement procedure of J.P. SELECTA, s.a.*

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INSTALLATION

Put the sterilizer on a flat, horizontal, level surface leaving a space of about 10cm at the back of the equipment.

For a better sterilizer function, do not situate it in places where it can be directly exposed to the sun radiation or heat sources.

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Do not use the incubator to dry or for thermal treatment of products which can give off vapours capable of making an explosive mixture.

Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

Do not use the equipment if it is not earthed.

If you change the plug, bear in mind the following

**Blue cable: Neutral.
Brown cable: Phase.
Yellow/green cable: Earth.**

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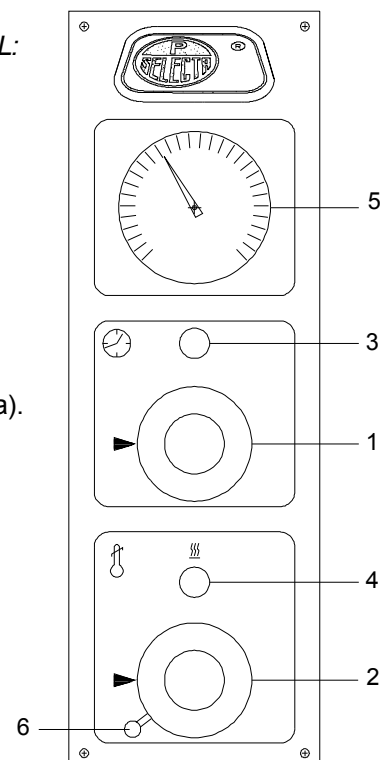
STARTING UP:

1. *Select the working temperature with the thermostat control knob (2).*
2. *Block the temperature control knob (2) with the blockage device (6) to avoid accidentally moving the knob.*
3. *Turn the main switch / switch selector (1) until the end of the scale. At the same time the mains indicator lamp (green) (3) and the heater operation indicator lamp (amber) (4) will light up.*
4. *Once the working temperature is reached, adjust the sterilization time turning the switch selector (1) at the desired time.*
5. *The thermostat scale is intended as a guide, the real temperature level inside the oven is the one that the thermometer shows.*

COMMENTS:

- * *Do not use the sterilizer at a lower temperature than ambient because the thermostat will not operate correctly.*
- * *All the temperature regulations must always be done in ascent. Before choosing any value, the knob must be turned left to the end of the scale. Then, a more exact regulation is obtained, as it compensates for the thermostat spindle inevitable movement.*

SUGGESTED STERILIZATION TIME WITH DRY HEAT:

140°C	2 hours
160°C	1 hour
170°C	40 minutes
180°C	20 minutes

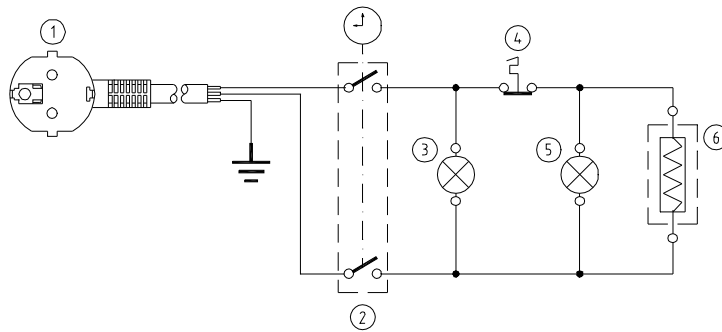
GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damages caused by an incorrect use or causes beyond the control of J.P. SELECTA, S.A.

Any manipulation of the equipment by personnel not authorized by J.P. SELECTA, S.A. automatically cancels the guarantee.

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MAINTENANCE

Before removing the casing, disconnect the equipment from the mains.

The manipulation of the internal electronic circuits of the incubator by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA, s.a. authorized technical services.

CLEANING:

For the cleaning of the different parts, use the following products:

Cleaning of stainless steel: alcohol.

Cleaning of plastic: alcohol with cotton duster.

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SPARE PARTS

To guarantee the safety of the equipment, all spare parts must be purchased from J.P.SELECTA, s.a.

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