**BOMBA MEMBRANA DE VACÍO  
DIAPHRAGM PUMP FOR VACUUM**

<b>BOMBA VACUM-SEL</b>	<b>3001001</b>
<b>VACUM-SEL PUMP</b>	<b>3001001</b>

**BOMBA MEMBRANA DE VACÍO Y PRESIÓN  
DIAPHRAGM PUMP FOR VACUUM AND PRESSURE**

<b>BOMBA VACUM-PRES</b>	<b>3001002</b>
<b>VACUM-PRES PUMP</b>	<b>3001002</b>

**INFORMACIÓN GENERAL****GENERAL INFORMATION**

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*The standard equipment consist of the following components:*

<b>Pieza</b>	<b>Piece</b>	<b>Código / Code</b>	<b>Cant. / Qty.</b>
Bomba VACUM-SEL	VACUM-SEL pump	3001001	1
Bomba VACUM-PRES	VACUM-PRES pump	3001002	1
Manual de instrucciones	Instruction Manual	80074	1

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*Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.*

<b>Código / Code</b>		<b>3001001</b>	<b>3001002</b>
<b>VACÍO / VACUUM</b>	<b>Caudal l/minuto / Flow rate l/minute</b>	1,08	
	<b>Vacío límite / Vacuum limit</b>	-0,6 bar ≈ 440mm Hg	
<b>PRESIÓN / PRESSURE</b>	<b>Caudal l/minuto / Flow rate l/minute</b>	--	1,3
	<b>Presión máxima / Maximum pressure</b>	--	1,2 bar
<b>Medidas exteriores cm / Overall dimensions cm</b>	<b>Alto / Height</b>	19	
	<b>Ancho / Length</b>	14	
	<b>Fondo / Width</b>	19	
<b>Consumo W / Consumption W</b>		40	
<b>Peso Kg / Weight Kg</b>		3	3,2

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*VACUM-PRES and VACUM-SEL pumps are equipped with a teflon member, housing and viton valves.*

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*It has a scale pressure gauge and a scale vacuum gauge according to the model and vacuum/pressure regulation valve (VACUM-PRES).*

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## INSTALLATION

Place the pump on a flat horizontal, level surface, leaving a space at the back and sides of the apparatus of about 10cm.

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Do not use the apparatus to dry or for thermal treatment of products which can give off vapours capable of making an explosive mixture.

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THE VACUM-SEL VACUM-PRES PUMPS ARE NOT SUITABLE FOR DECATING (MIXING) LIQUIDS.

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A VACUUM TRAP IS HIGHLY RECOMMENDABLE IN THE CONNECTION TO THE PUMP TO AVOID THE ACCIDENTAL LEAK OF LIQUIDS.

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Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

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Do not use the apparatus if it is not earthed.

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If you change the plug bear in mind the following:

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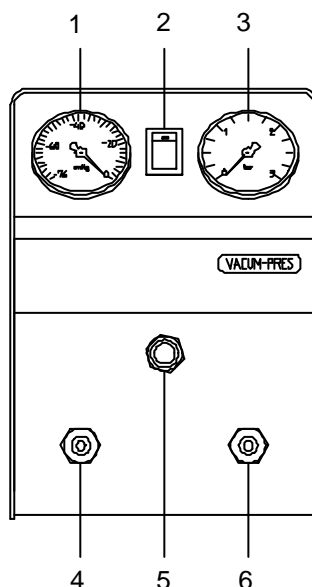
Blue cable: Neutral.  
Brown cable: Phase.  
Yellow/green cable: Earth.

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
CONTROL PANEL:

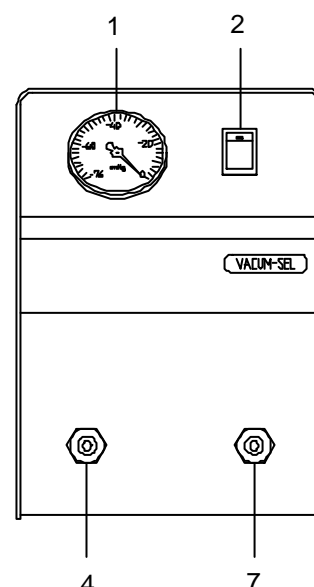1. Scale vacuum gauge from 0 to -1 bar.
2. Main switch.
3. Scale pressure gauge from 0 to 2.5 bar.
4. Vacuum inlet.
5. Pressure / vacuum regulator.
6. Pressure outlet.
7. Air outlet.

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**STARTING UP:**

1. Connect the tubes to the inlet / outlet 4, 6 and 7.
2. Press the main switch (2).
3. Control the pressure / vacuum with the scale pressure gauge (1) / scale vacuum gauge (3).
4. Regulate the pressure / vacuum desired with the regulator (5).

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**MAINTENANCE**

Before removing the cover disconnect the apparatus from the mains.

The manipulation inside of the equipment by unauthorized personnel can cause irreparable damage. Take it to one of the J.P. SELECTA authorized technical services.

**CLEANING:**

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with cotton duster.

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**SPARE PARTS**

To guarantee the safety of the equipment, the spare parts must be purchased from J.P. SELECTA, S.A.

Descripción	Description	3001001	3001002
1. Tetina toma de vacío	Vacuum inlet	15043	15043
2. Tetina expulsión aire	Air outlet	15050	-
3. Tetina toma presión	Pressure inlet	15052	15052
4. Vacuómetro	Scale vacuum gauge	16012	16012
5. Manómetro	Scale pressure gauge	-	16011
6. Interruptor	Switch	20034	20034
7. Microbomba	Pump	24204	24204
8. Regulador de flujo	Flux regulator	-	34040

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## GUARANTEE

*This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA, S.A.*

*Any manipulation of the apparatus by personnel not authorized by J.P. SELECTA, S.A. automatically cancels the guarantee.*

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