U.V. ESTÉRIL

3000416

**PUPITRE DE ESTERILIZACIÓN POR RAYOS ULTRAVIOLETA
ULTRAVIOLET RADIATION STERILIZATION DESK**

INFORMACIÓN GENERAL

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without first reading this handbook.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT!**
J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.
- 6) *If any modification, elimination or lack of maintenance of any device of the equipment by the user transgresses the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

INSTALLATION

Before connecting the apparatus into mains, make sure that the voltage indicated on the plate at the back of the apparatus corresponds to the mains voltage.

On the frontal part is placed the main switch with indicator lamp which starts up the ultraviolet lamp.

Due to the intense short wave ultraviolet radiation, take care to protect eyes and skin from direct radiation. For this purpose, the apparatus is provided with a microswitch which disconnects the lamp on opening the sterilizer cover.

The ultraviolet rays are radiated by a germicidal lamp of 15W or 8W depending on model and with a wavelength of 2.537 Angstroms.

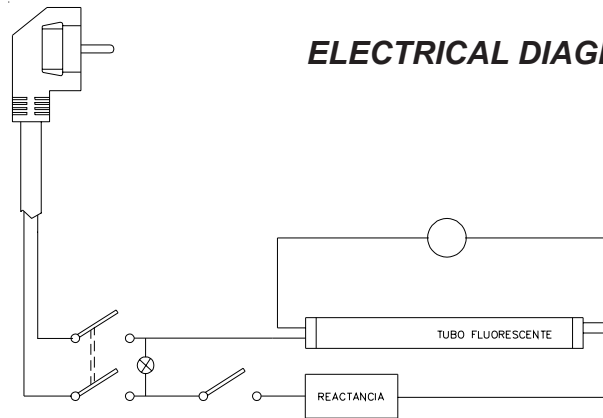
STERILIZATION TIME:

For clinical and odontological instruments, Petri capsules, Staphylococcus, Streptococcus, Bacillotific, contaminations, etc. five minutes radiation is sufficient. For B. Subtilis, Penicium Glaucum, Phlei resistant acid, spores type contamination sterilization is obtained in 15 minutes of radiation.

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The life of an ultraviolet bulb is approximately 5,000 hours.

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IMPORTANT

This apparatus should in no way substitute the classical means of sterilization such as autoclaves, poupinel, etc... It is normally used to maintain previously sterilized materials for seeding, etc... Anyway, germicidal lamps can eliminate some kind of bacteriums, moulds and algae.

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GUARANTEE

This product is guaranteed for one year. The guarantee does not cover damage caused by incorrect use or causes beyond the control of J.P. SELECTA, S.A.

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Any manipulation of the apparatus by personnel not authorized by J.P. SELECTA, S.A. automatically cancels the guarantee.

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*Declares that the equipment:
Model: U.V ESTERIL Code: 3000416*

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*73/23/CEE Low tension regulation
89/336/CEE Electromagnetic compatibility regulations*

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By fulfilling the following International Standards:

EN 50081-1 EN 50082-1 EN 61010-1

EN 50081-1 EN 50082-1 EN 61010-1