PLACA CIRCULAR	1000718
PLACA CIRCULAR	1000442
PLACA CIRCULAR	1000443
SELECTEMP	1000444
PLACFIN	7000497
UNIPLAC	6002312
UNIPLAC	6002314
UNIPLAC	6002316

**PLACAS CALEFACTORAS
HOTPLATES**

INFORMACIÓN GENERAL

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GENERAL INFORMATION

- 1) *Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.*
- 2) *Do not install or use the equipment without reading this handbook before.*
- 3) *This handbook must always be attached to the equipment and it must be available for all users.*
- 4) *If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta's technical service.*
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) *If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.*
- 7) *Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.*

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CONTENTS

GENERAL INFORMATION
CONTENTS
PACKING LISTS
ACCESSORIES
EQUIPMENT DESCRIPTION
TECHNICAL FEATURES
INSTALLATION
OPERATION
MAINTENANCE
SPARE PARTS
ELECTRICAL DIAGRAM
GUARANTEE
"EC" CONFORMITY DECLARATION

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The standard equipment consist of the following components:

Descripción	Description	Código / Code	Cant. / Qty.
Placa circular	<i>Circular hotplate</i>	1000718	1
Placa circular	<i>Circular hotplate</i>	1000442	1
Placa circular	<i>Circular hotplate</i>	1000443	1
Placa circular "Selectemp"	<i>Circular hotplate "Selectemp"</i>	1000444	1
Placa rectangular "Placfin"	<i>Rectangular hotplate "Placfin"</i>	7000497	1
Placa circular "Uniplac"	<i>Circular hotplate units "Uniplac"</i>	6002312	1
Varillas verticales Ø12x60	<i>Vertical rods Ø12x60</i>	12103	2
Varilla transversal Ø12x45	<i>Horizontal rods Ø12x45</i>	12105	1
Nuez sujeta varillas	<i>Bosshead for rods</i>	7008133	2
Llave allen	<i>Head key</i>	9100005	1
Placa circular "Uniplac"	<i>Circular hotplate units "Uniplac"</i>	6002314	1
Varillas verticales Ø12x60	<i>Vertical rods Ø12x60</i>	12103	2
Varilla transversal Ø12x77	<i>Horizontal rods Ø12x77</i>	12101	1
Nuez sujeta varillas	<i>Bosshead for rods</i>	7008133	2
Llave allen	<i>Head key</i>	9100005	1
Placa circular "Uniplac"	<i>Circular hotplate units "Uniplac"</i>	6002316	1
Varillas verticales Ø12x60	<i>Vertical rods Ø12x60</i>	12103	2
Varilla transversal Ø12x105	<i>Horizontal rods Ø12x105</i>	12107	1
Nuez sujeta varillas	<i>Bosshead for rods</i>	7008133	2
Llave allen	<i>Head key</i>	9100005	1
Manual de instrucciones	<i>Instruction manual</i>	80061	1

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EQUIPMENT DESCRIPTION

The hotplates are designed to work with a big temperature range, up to 130°C (Placfin), 250°C (Selectemp) up to 400°C (Circular hotplates and Uniplac).

Temperature regulation is done by a hydraulic thermostat (Selectemp and Placfin) or impulse regulators (Uniplac) independent for each plate.

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TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

		Placas circulares / Circular hotplates			Placa circular / Circular hotplate SELECTEMP	Placa rectangular / Rectangular hotplate PLACFIN	Placas circulares / Circular hotplate units UNIPLAC		
Código / Code		1000718	1000442	1000443	1000444	7000497	6002312	6002314	6002316
Nº de plazas / Nº of plates		1	1	1	1	1	2	4	6
Medidas útiles de la placa cm. / Plate dimensions cm	Ancho / Length	--	--	--	--	23	--	--	--
	Fondo / Width	--	--	--	--	22	--	--	--
Ø Placa en cm. / Ø Plate cm		12	15	19	15	--	12		
Medidas exteriores cm. / Overall dimensions cm	Alto / Height	8			6	8	23		
	Ancho / Length	20	22	26	30	22	47	77	105
	Fondo / Width	20	22	26	22	30	27		
Consumo W. / Consumption W		1000	1500	2000	1500	400	1000	2000	3000
Temperatura máxima / Maximum temperature		400°C			250°C	130°C	400°C		
Peso Kg / Weight Kg		2.1	3	4.5	3	3.7	8	15	22

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INSTALLATION

Place the hotplate on a flat horizontal, level surface, leaving a space at the back and sides of the apparatus of about 10cm.

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For a longer life of the apparatus provided with hotplate, it is recommended that its surface be covered as much as possible by the beaker to be heated. This avoids overheating and deformation of the plate by absorbing more the heat.

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Because of that, when using round-bottomed, Soxhlet or Kjeldahl flasks it is necessary to use the adapters in aluminium described in "HOTPLATES" section of our catalog.

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Be sure that the voltage supply is the same as the one indicated on the characteristics plate.

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Do not use the apparatus if it is not earthed.

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If you change the plug bear in mind the following:

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*Blue cable: Neutral
 Brown cable: Phase
 Yellow/green cable: Earth*

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
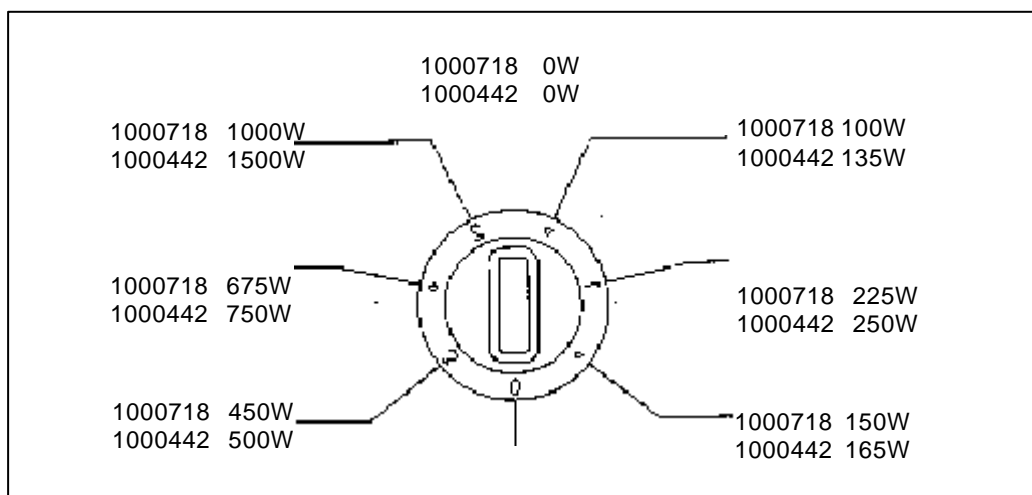
These heaters include a selector of 6 heating powers settings and stop. In the table below the power in watts is given according to the selector position.

PANEL DE MANDOS PLACAS CIRCULARES EN BATERIA UNIPLAC:

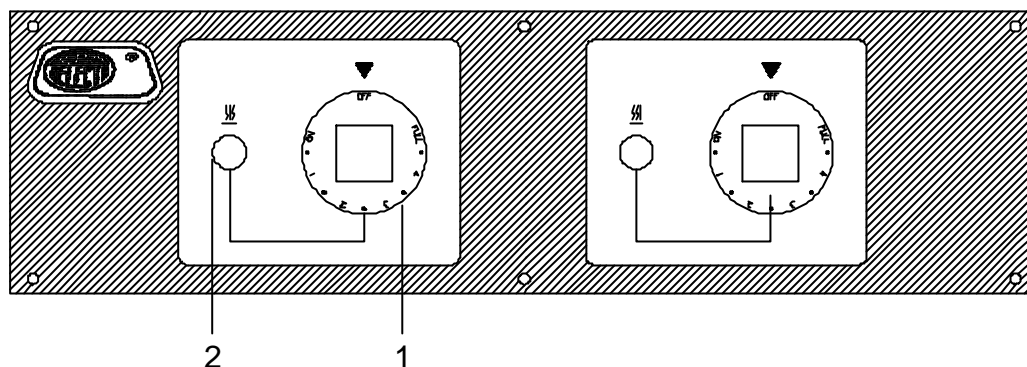CIRCULAR HOTPLATE UNIPLAC UNITS CONTROL PANEL:

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CIRCULAR HOTPLATE UNIPLAC UNITS STARTING UP:

1. Move the control knob of temperature (1) from the "OFF" position (disconnected), to the "FULL" position (maximum power).
2. It will be noted that according to the control position, the ON-OFF cycle switch of the indicator lamp for the working of the heater (2) will be longer or shorter, which will permit adjusting to the desired working temperature.
3. To stop, move the control button to "OFF" position.

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CIRCULAR HOTPLATE SELECTEMP:

Circular hotplate "SELECTEMP" is provided with a temperature regulator thermostat whose sensitive part is in the center of the plate in order to register the temperature of the container to be heated. Move the thermostat button numbered 1 to 12 to the right and fix it at one point, bearing in mind that the higher the point fixed, the higher the temperature. If after a reasonable time the desired temperature has not been reached, move the button of the thermostat to more or less, as desired. To disconnect turn the button to 0.

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IMPORTANT

Treat the central thermostat of the plate very carefully, do NOT try to scrape it with anything to clean it. Should any liquid spill onto the plate simply clean it with a damp cloth.

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Follow the same instructions as for "SELECTEMP" plate and read the temperature with the thermometer included in the apparatus.

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WARNING!!! Caution with the hotplate surface, serious burns can be produced, even after switching off the apparatus.

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MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

The manipulation of the internal electronic circuits of the apparatus by unauthorized personnel can cause irreparable damage. Take it to one of the J.P.SELECTA authorized technical services.

CLEANING:

For the cleaning of the different parts use the following products:

Cleaning of stainless steel: Alcohol

Cleaning of plastic: Alcohol with cotton duster.

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Any manipulation of the apparatus by unauthorized personnel by J.P. SELECTA, S.A. cancels the guarantee automatically.