## REACTOR PARA PLANTA PILOTO REACTOR FOR PILOT PLANT

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# INDEX

<b>1. General .....</b>	36
1.1 Introduction .....	36
1.2 Safety.....	36
1.3 Equipment Description.....	39
1.4 Equipment Configuration.....	40
1.5 Control panel.....	41
<b>2. Unpacking and Installation .....</b>	42
2.1 Prepare .....	42
2.2 Installation .....	42
2.3 Electric control system connection .....	49
2.4 Connect the glass reactor with vacuum pump and temperature control unit .....	49
2.5 Operation .....	54
<b>3. Technical Data .....</b>	55
3.1 Borosilicate Glassware 3.3 .....	55
3.2 Lubrication.....	56
<b>4. Operating .....</b>	57
4.1 Caution .....	57
4.2 Operating .....	58
<b>5. Maintenance and Management.....</b>	59
5.1 Management.....	59
5.2 Cleaning.....	59
5.3 Maintenance .....	59
5.4 Troubleshooting .....	60
<b>6. Accesories and Spare Parts.....</b>	61
<b>7. Examples of Use of Support Trolleys .....</b>	66
<b>8. Application examples .....</b>	67

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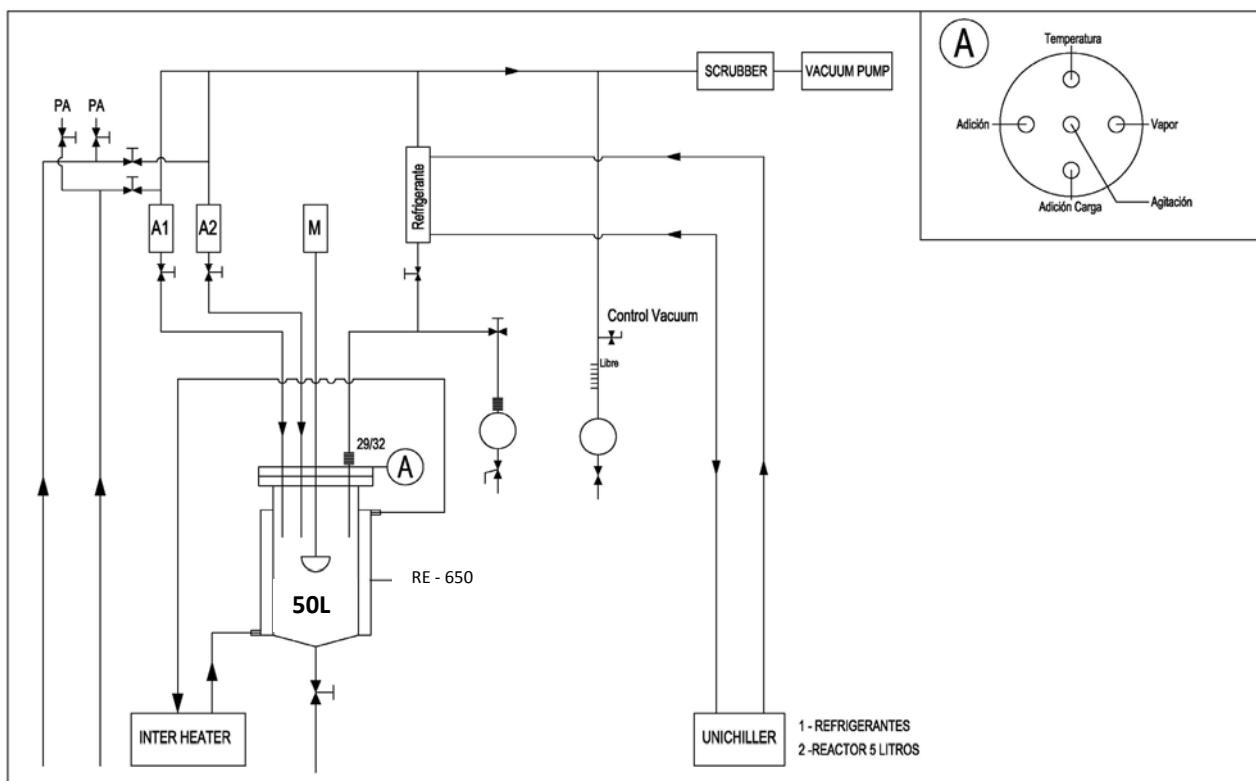
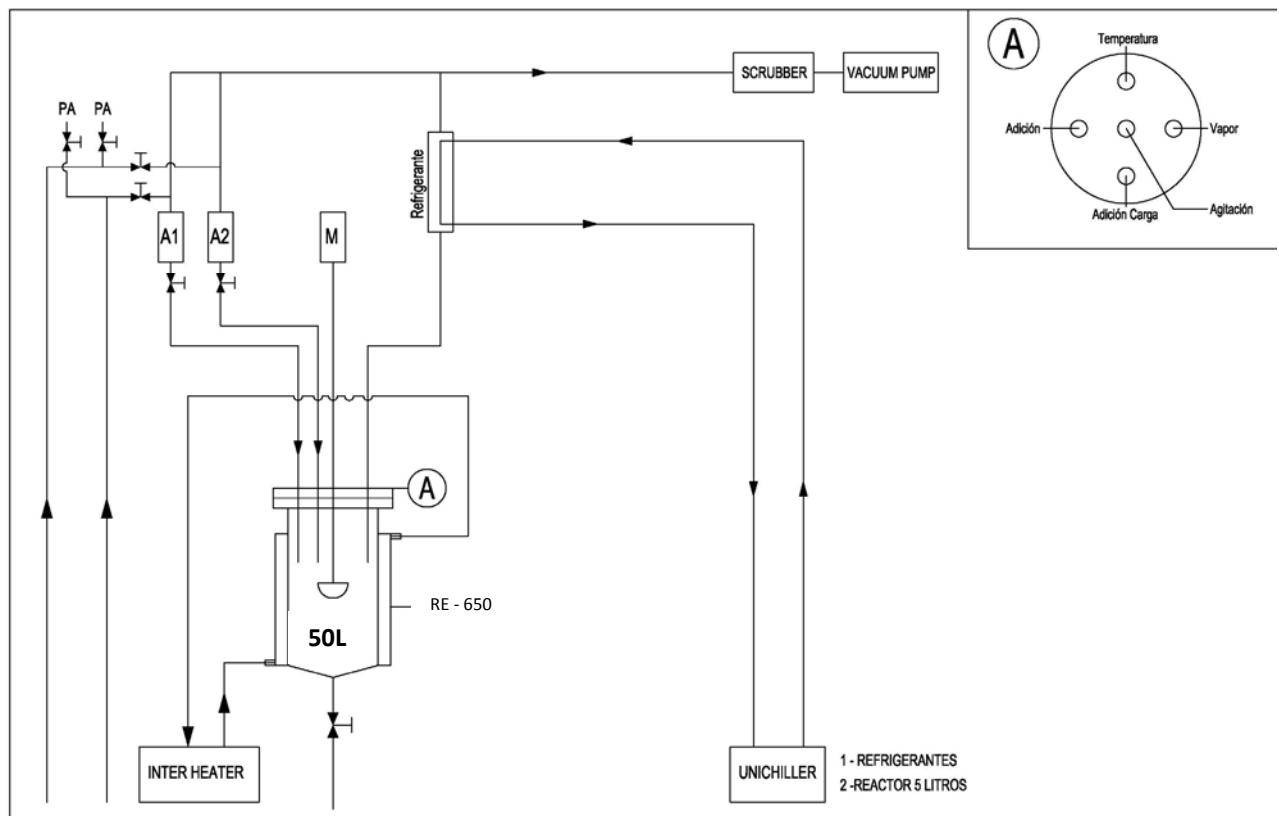
# **1. General**

## **1.1 Introduction**

Thanks for using 'GR-series speed-regulation glass reactor'. This operation manual stated the installation, using, maintaining and servicing of glass reactor. Before using this instrument, please ensure that you have read and understood this manual completely.

## **1.2 Safety**

This chapter describes the installation, safety rules in the process of using of 'GR-series speed-regulation glass reactor'. Users must grasp the related warning signs, strictly abide by the operation procedures to ensure the security of the equipment and personal and avoid the occurrence of accident.

### **1.2.1 User's Qualification**

GR-series speed-regulation glass reactor must be operated by the person who has the practical operating experiences and can grasp of the detailed requirements in this manual. Otherwise, it must be used under the guidance of the person who has the related technology skills.

### **1.2.2 Proper Use**

Application fields:

- a) The process which need high temperature, low temperature conditions.
- b) The process which need reduce pressure and stirring at the same time.
- c) Dispersion process.
- d) Distillation (reflux), concentration.
- e) Assembled into a glass reactor total system according to the specific requirements.

### **1.2.3 Improper Use**

The operation that not abide by the related stipulation in this manual are regarded as improper use. Any damage caused by improper use is responsible by the users themselves.

Operation under the following conditions is prohibited:

- a) Used in explosive gas environment or explosive dust environment.
- b) Used in the places where the power supply is not in conformity with the requirements.
- c) Used in high magnetic fields, corrosive environment.
- d) Hard, brittle bulk materials (such as stones, soil samples, etc.) may damage the vessel.
- e) The volume of the sample added to the vessel exceeds the specified limit.

#### 1.2.4 Warning Sign

Sign	Description
	<b>Danger</b> It shows that the situation is very dangerous and will lead to death or serious injury.
	<b>Warning</b> It shows that the situation is very dangerous and will lead to death or serious injury.
	<b>Prohibit</b> Get wet in the rain or splashing water.
	<b>Note</b> It shows that the situation is very dangerous and will lead to injury.
	<b>Note</b> Caution high temperature.
	<b>Note</b> Please wear protective equipment, otherwise may cause personal injury.
Reminder	<b>Reminder</b> May cause equipment damage.

Especially pay attention to each warning sign.

### 1.2.5 Hazards Related to the Instrument

Please pay attention to the following safety tips:

 Danger	<b>Danger</b> <ul style="list-style-type: none"><li>● Do not use it in explosive gas environment or explosive dust environment.</li></ul>
 Warning	<b>Warning</b> <ul style="list-style-type: none"><li>● Make sure the power supply is in conformity with the requirements on the name plate</li><li>● Equipotential connection has been done, do not remove inner grounding connection line.</li><li>● Before open the equipment enclosure for maintenance and repair, be sure to disconnect the power supply, then operate it after 5s, ensure the residual voltage release to safety value to avoid electric shock.</li><li>● Do not use broken glass parts</li></ul>
 Prohibit	<b>Prohibited</b> <ul style="list-style-type: none"><li>● Do not use it outdoor. Get wet in the rain or splashing water will cause electrification on metallic shell surfaces, and will cause casualties.</li></ul>
 Warning	<b>Warning</b> <ul style="list-style-type: none"><li>● Do not touch the vessel surface and connection pipes with your body directly to avoid scalding.</li></ul>
 Note	<b>Note</b> <ul style="list-style-type: none"><li>● If there's any abnormal situation when using it, please disconnect the power supply immediately for troubleshooting or contact professional maintenance personnel.</li><li>● Be aware of broken glass parts.</li><li>● It may have static electricity when adding material to the vessel through feeding port</li></ul>
Reminder	<b>Reminder</b> <ul style="list-style-type: none"><li>● Do not use it in the environment mentioned above. Otherwise, it will affect its normal running and lifetime.</li><li>● Keep well ventilated around the equipment.</li><li>● Do not use it when there's no one there.</li></ul>

### 1.2.6 Other Hazards

 Warning	<ul style="list-style-type: none"><li>● The sample in the vessel or some solvents near it may form peroxides, high concentrations of flammable solvents, etc.</li><li>● Acid, alkali steam and other corrosive materials around the equipment will damage the equipment insulation and affecting component performance and service life.</li><li>● Beware of the risk of explosion when dealing with hazardous materials or unknown samples.</li></ul>
--	--

### 1.2.7 Safety Measures

 	<ul style="list-style-type: none"><li>● Please wear personal protective equipment when operating this instrument, such as protective glasses, protective clothing and gloves</li></ul>
---	--

---

### 1.3 Equipment Description

RE-650 speed-regulation glass reactors are mainly used for synthetic reaction, distillation and concentration of different types of materials. We have single wall, double wall and triple wall glass reactors.

The reactor can be pumped to a negative pressure state according to some experimental requirements. A constant pressure funnel or an adjusting valve on the feeding bottle allows you to add material to the vessel at a uniform and controlled speed. Condenser helps to recover some materials distilled during reaction.

Single wall glass reactor has no jacket.

Double wall glass reactor has a jacket. Circulating liquid can be filled into the jacket to heat or cool materials inside.

Triple wall glass reactor has two jackets, the inner jacket is for circulating liquid, and the outer layer should be pumped to vacuum condition for thermal insulation.

## 1.4 Equipment Configuration

RE-650 speed-regulation glass reactor (Customized product as per the real object)

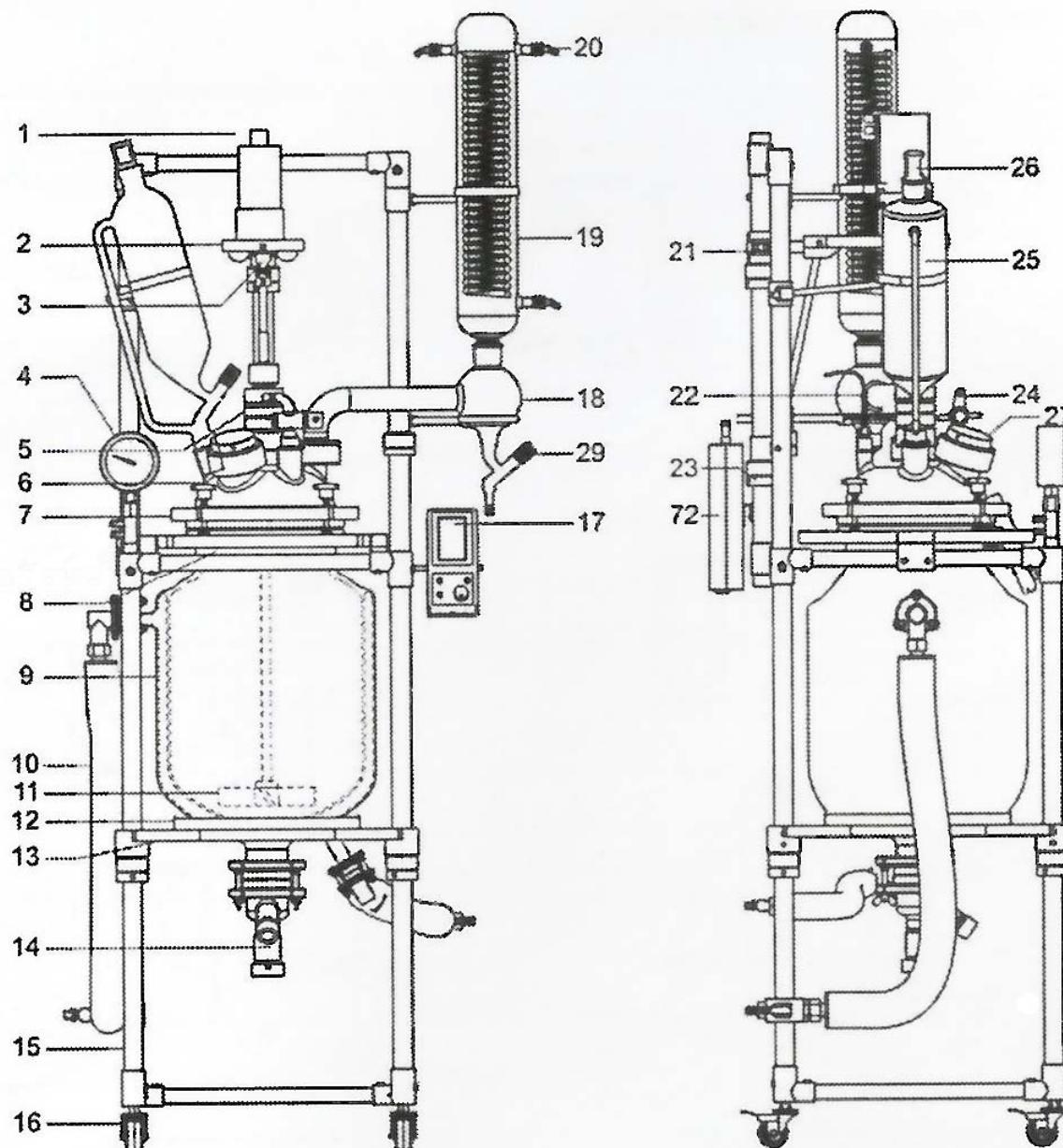
Figure 1.4.2 - RE-650 Speed-regulation glass reactor

1 Motor	2 Motor support	3 Universal coupling	4 Vacuum gauge
5 Stirrer bearing guide	6 Reactor lid	7 Upper sealing clamp	8 Lower sealing clamp
9 Reactor vessel	10 Circulating hose	11 Stirring paddle	12 Tray gasket
13 Vessel supporting platform	14 Discharge valve		
15 Frame	16 Caster	17 Control box	18 Condensing dispenser
19 Condenser	20 Joint	21 Fixing frame	22 Temperature sensor
23 Fine tuning parts	24 Feeding valve	25 Constant-pressure funnel	26 Plug
27 Charging port sealing kit	29 15# cock	72 Motor controller	


## 1.5 Control panel

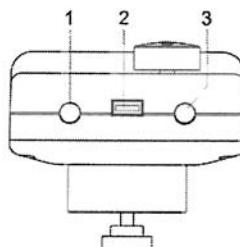Control panel

1 Mode	2 Set
3 Power on/off	4 Start/stop
5 Stirring time	6 Temperature
7 Torque	8 Fault information
9 Actual stirring speed	10 Setting stirring speed


5. MODE: monitoring interface. Press this button to view fault information. Press this button again to go back to monitoring interface. Press the button longer to enter password setting interface (parameters have been set ex-factory).
6. SET: Monitoring interface. Press this knob to enter stirring speed setting mode. Turn the knob slowly to target value and press it again to save the setting speed. If you do not press the button again, the setting speed will not be saved.
7. POWER ON/OFF: Power on and off
8. START/STOP: Start or stop stirring

Interfaces on the control panel

1 Temperature sensor socket    2 USB interface    3 485communication interface (optional)

**Motor controller**

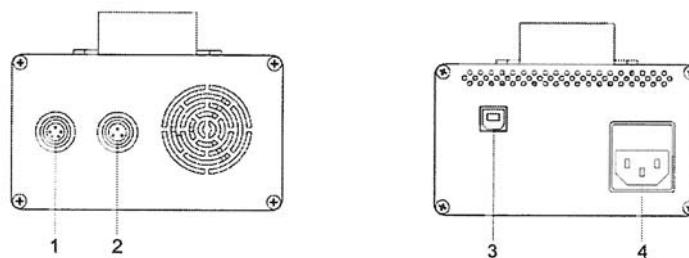Interfaces on the motor controller

1 Signal interface	2 Motor power supply socket
3 USB Interface	4 Main power supply socket

## 2. Unpacking and Installation

### 2.1 Prepare

- 1) After unpacking the case, please read the operation manual carefully, and check whether the components are enough compare with the packing list.
- 2) Take out the frame, put it on the floor then lock the casters.

Note: The vessel and frame are assembled completely of reactors less than 20L.

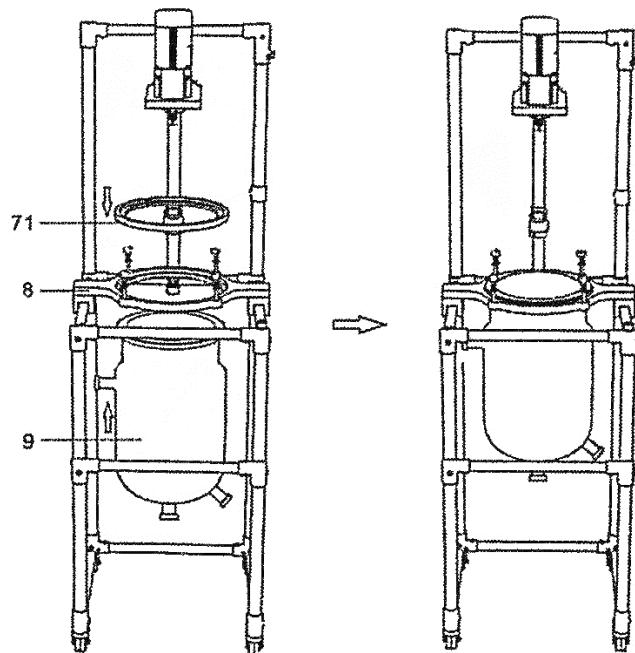
- 3) Please clean the glass parts to maintain its cleanliness to meet the test requirements before installation.

Note: Glass device is fragile, please be careful when unpacking.

### 2.2 Installation

#### 2.2.1 RE-650 vessel installation

- 1) The vessel (9) fits from the bottom of the middle tray (8).
- 2) The sealing ring (71) is fitted from the flange mouth of the vessel (9) .
- 3) Put the sealing ring (71) on the middle tray (8) to make their circle fitted.

**Reactor Vessel Installation**

8 Lower sealing clamp

9 Reactor vessel

71 Lid gasket

## 2.2.2 Stirring paddle installation

Four plum knobs loose to the bolts can be flat, remove the upper tray (7), tray gasket, reactor lid (6) and reactor lid openings sealing gasket. Place the stirring paddle (11) in the vessel body. Coated vacuum grease on the contact surface of the vessel body and the reactor lid. Install the reactor openings sealing gasket, reactor lid (6), tray gasket (7) and upper tray (7) in turn, tighten the plum knob.

Note: Please refer to 1.4 for the mounting position of the reactor lid.

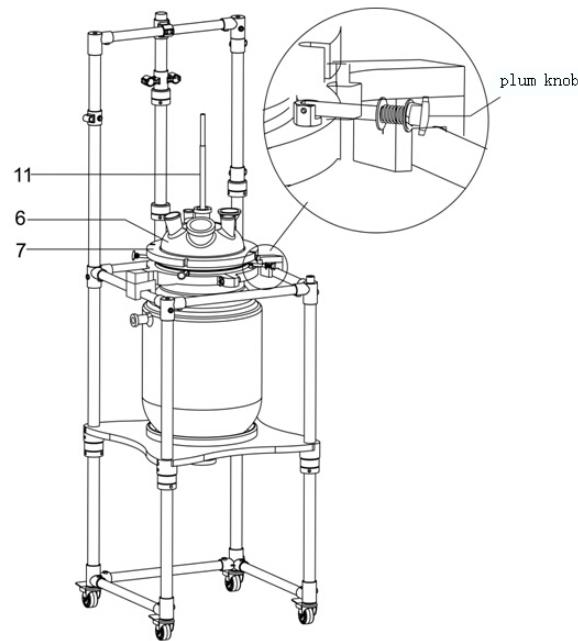
Figure 2.2.2 Stirring paddle installation

6 Reactor lid      7 Upper tray      11 Stirring paddle

### 2.2.3 Stirring shaft installation

Please install the stirring paddle according to figure 2.2.3.

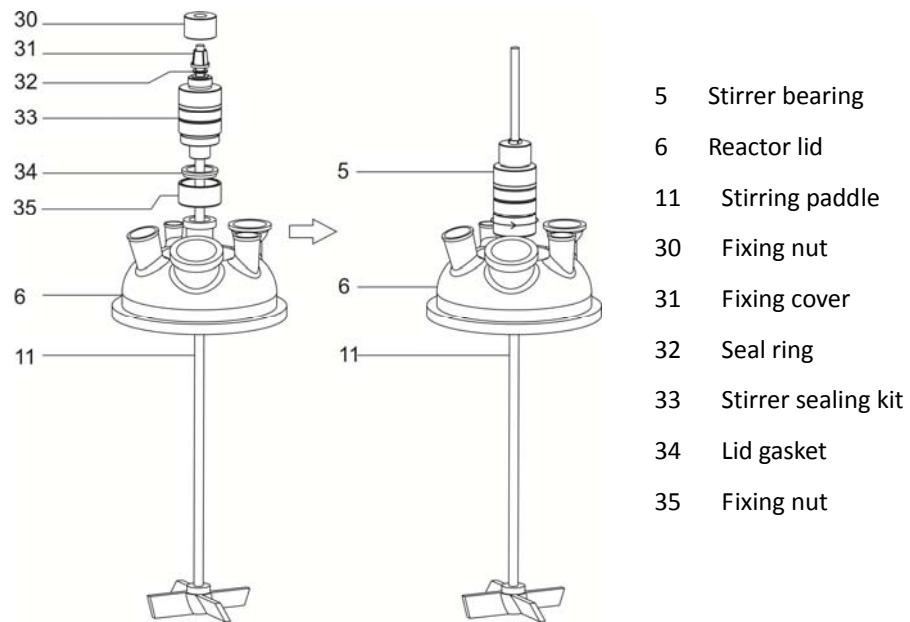
Figure 2.2.3 Stirrer bearing installation

### 2.2.4 Motor installation

The motor, motor tray, universal coupling and fixing pole have been installed together before leaving factory, please see figure 2.2.4.1.

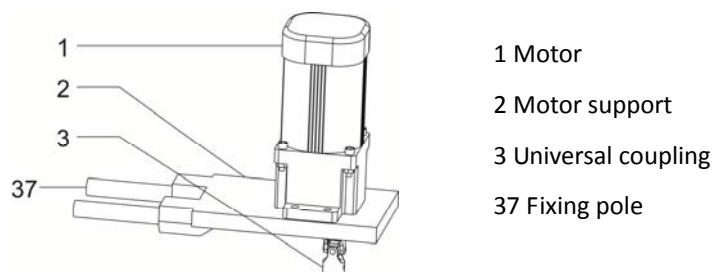
Figure 2.2.4.1 Motor assembly

---

### 1) RE-650 motor kit installation

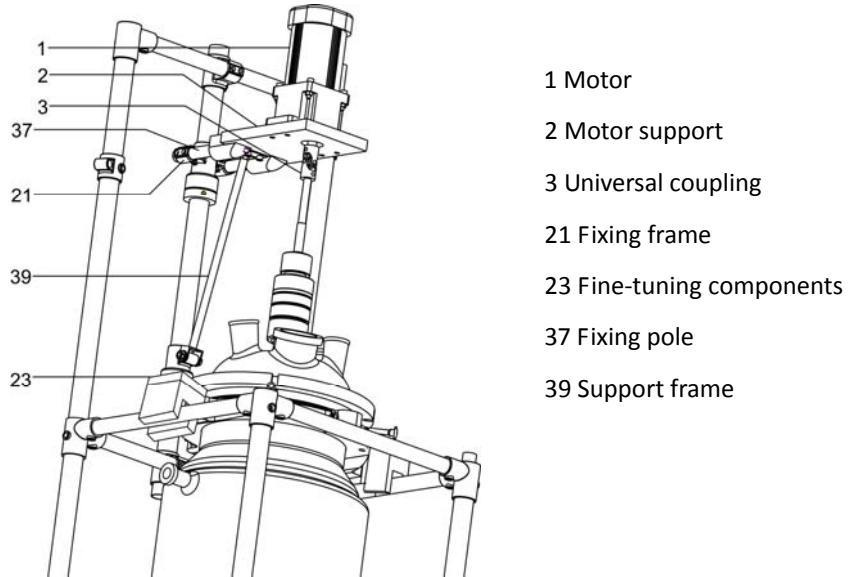
Figure 2.2.4.3 RE-650 motor installation

Connect support frame (39) and motor support (2), tighten the connection screw. Install the motor kit on the fixing frame (21) to make the motor axis and stirring paddle coaxial, fix the universal coupling (3). Adjust the position of fixing frame (21) to make the gap between bottom of the stirring paddle and vessel bottom >30mm, fix the fixing frame (21). Then tighten the connecting bolt and screw on the lower position of the support frame (39). Adjust the fine-tuning components (23) to make the level of the motor tray is horizontal. Tighten the fixing nut (30), fix the stirring paddle.

### 2.2.5 Installation of circulation connectors of the condenser

Please install it according to figure 2.2.5.

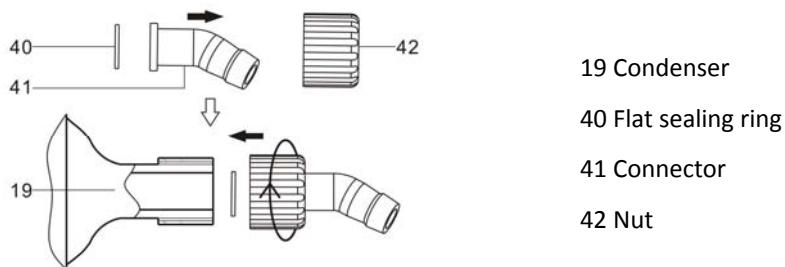
Figure 2.2.5 Nozzle and condenser installation

## 2.2.6 Condensing dispenser and condenser installation

Fix the support (44) on the frame, put the condensing dispenser (18) on the support (44). Connect the condensing dispenser (18) and the reactor lid joint, adjust the position of support (44), tighten the fixing screw (45) and fixing nut (48).

Put the condenser (19) on the upper joint of the condensing dispenser (18), the rubber bracket (43) is locked around the upper portion of the condenser (19), please see figure 2.2.6.2.

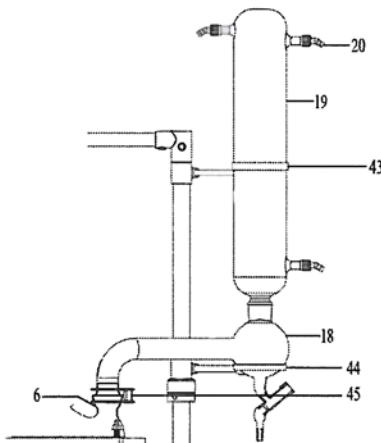
Figure 2.2.6 Installation of Condensing Kit

6 Reactor lid	18 Condensing dispenser	19 Condenser	20 Connector
43 Rubber bracket	44 Condensing dispenser support	45 Clamp	

## 2.2.7 Temperature sensor installation

Put the fixing screw (49), fixing cover (50), O-ring (51) and fixing base (70) on the sensor in turn, adjust sensor (22) position, tighten fixing nut (49).

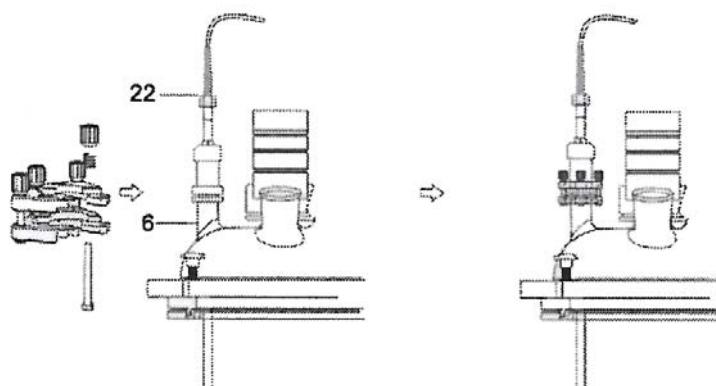
Figure 2.2.7 Installation of Temperature Sensor

6 Reactor lid	22 Temperature sensor
---------------	-----------------------

## 2.2.8 Solid charging port sealing kit installation

Please install it according to figure 2.2.8.

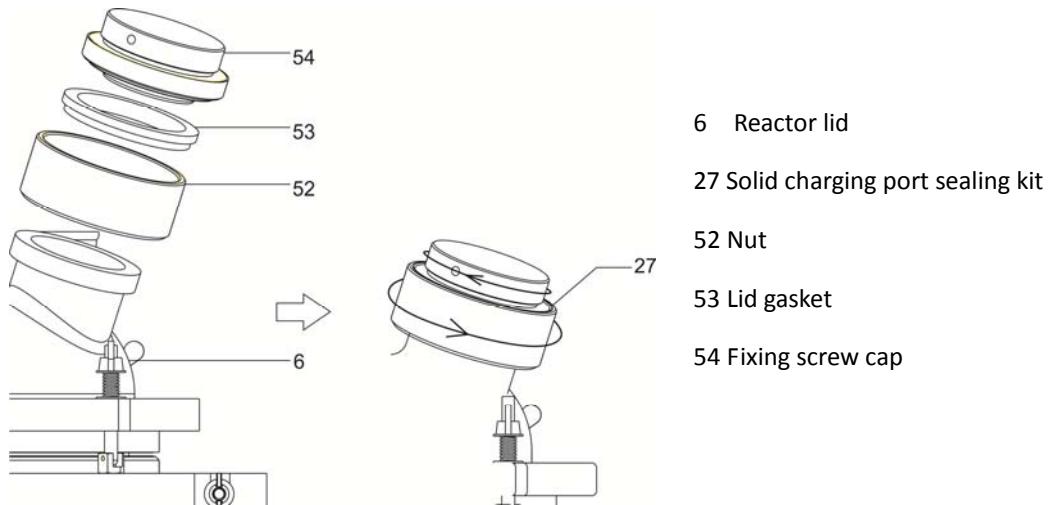
Figure 2.2.8 Solid charging port sealing kit installation

## 2.2.9 Constant-pressure funnel and feeding port valve installation

Please install it according to figure 2.2.9.

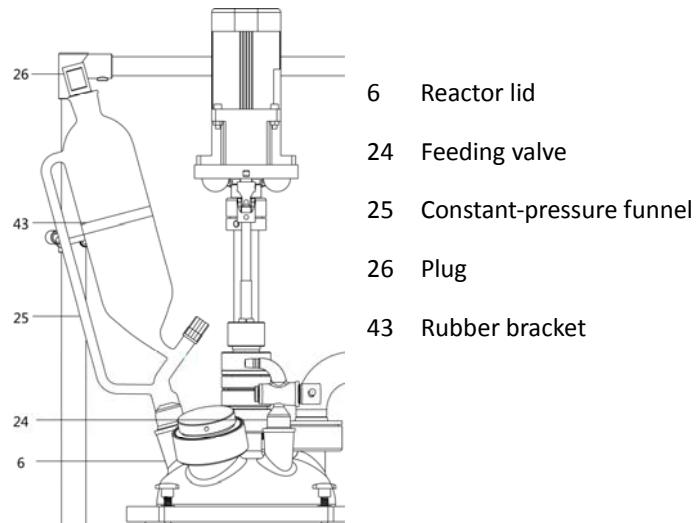
Figure 2.2.9 Constant-pressure funnel and feeding valve installation

## 2.2.10 Discharge valve installation

Note: Make sure that the discharge valve is open before installation, in case of damage of the reactor body.

Please refer to Figure 2.2.10.

Fit upper part of the flange (59) and a lid gasket (60) to the glass drain port (9) and the lower part of the flange and gasket to the discharge valve (14). Put sealing ring (61) between the upper and lower flange, keep them in the right position and tighten the butterfly nut.

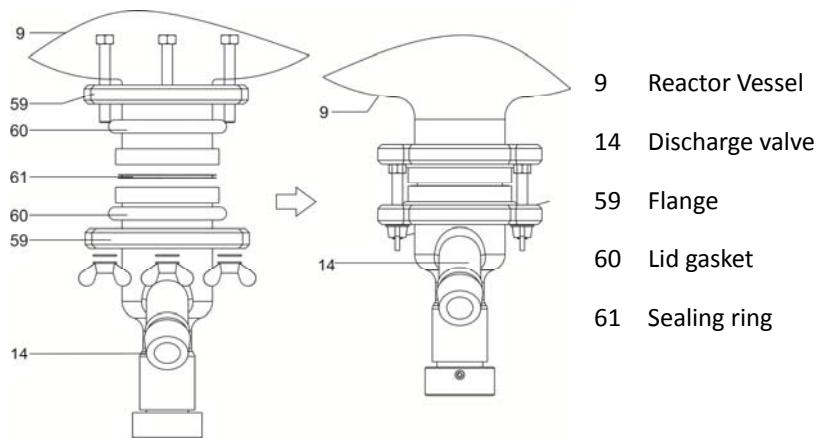
Figure 2.2.10 RE-650 discharge valve installation

## 2.2.11 Vacuum gauge and control box installation

Please see chapter 1.4 to find the installation position of Vacuum gauge (4) and control box (17).

## 2.2.12 Vessel jacket circulating pipe kit installation

The lower connecting joint of the vessel jacket is inlet, the upper is outlet. Please refer to figure 2.2.10.2 to install it.

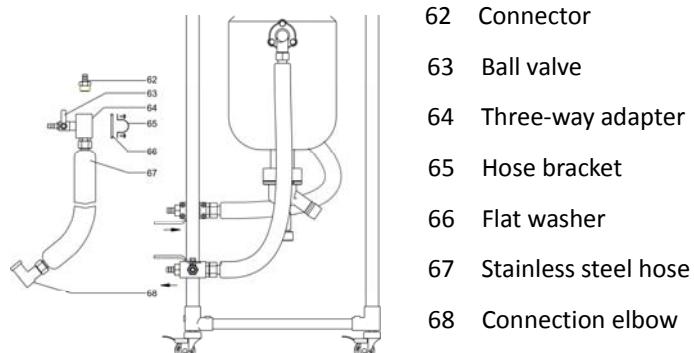
Figure 2.2.12 Circulating pipe kit installation

## 2.3 Electric control system connection

Connect the motor, the motor controller and the control box.

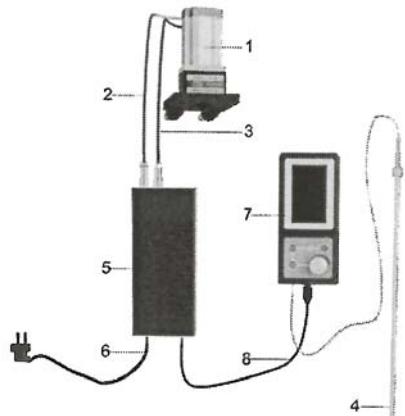
Figure 2.3 Connection of Moto /Motorcontroller/Control box

1 Motor	2 Signal detection wire	3 Motor power supply cord
4 Temperature sensor	5 Motor controller	6 Power supply cord
7 Control box	8 USB connection line	

## 2.4 Connect the glass reactor with vacuum pump and temperature control unit

Connect the glass reactor with Dynamic Temperature Control System and vacuum pump, make sure that there's no leakage between each joint.

Note: There are two ports on the back of the vacuum gauge. Connect one of the ports to the vacuum port of the condenser and the other to the vacuum pump.

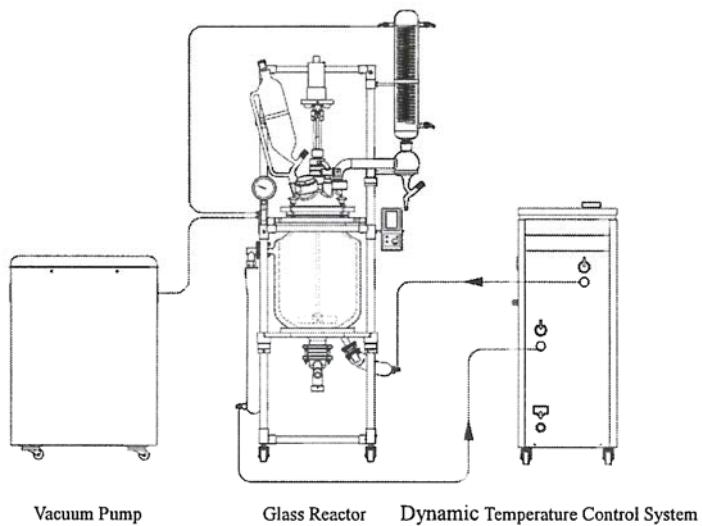
Figure 2.4 Diagram of Circulation Hose Connection

## Rod stirrers

FOR MIXTURES, HOMOGENIZATIONS, SUSPENSIONS, GASIFICATION OF SOLUTIONS AND RECIRCULATION OF SUBSTANCES.  
DESIGNED FOR CONTINUOUS WORK.

BRUSHLESS MOTORS FREE OF MAINTENANCE AND SPARKS.

ENCAPSULATED CASES PARTIALLY SEALED AGAINST AGGRESSIVE SUBSTANCES AND VAPOURS. THERMAL PROTECTION FOR OVERHEATING WITH SELF-DISCONNECTION AND INDICATOR LED. POWER SUPPLY VOLTAGE: 110 / 220V - 50-60 Hz.  
HEIGHT ADJUSTABLE AGITATION ROD (see wide selection of rods and accessories).

## Hei – TORQUE stirrers

**Powerful**  
Stirring

Newest motor generation for maximum power at **minimum noise level**  
**- below 50 db**

The **intuitive touch-panel** made of glass for easy operation

Sealed housing, which complies with the high protection class IP 54, guarantees longevity and maintenance-free **24-hour operation** in an aggressive environment

An overtemperature sensor prevents **heat-up situations** particularly valuable in unattended continuous operation

Increased safety with individual performance control:  
Set the **intensity of the starting operation**, the maximum **rotational speed**, and the maximum **torque limit** according to your application

The sealed panel made of glass **increases the tightness** of the entire housing

**Safe start and stop of operation** via slide touch panel to avoid unintended stirring

Control panel  
Hei-TORQUE Value

Control panel  
Hei-TORQUE Precision

Coolant double effect DIN 29/32 L: 300mm

Addition funnel 250ml, compensated pressure DIN29 / 32

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## 2.5 Operation

---

1. Turn the power switch on and the temperature and the speed will be displayed.
  2. The speed is adjusted to 0, click on the Rotation key, slowly adjust the speed, the rotation of the agitation blade, and observe if any anomaly occurs.
  3. Pour the materials into the container through solid feed or other inputs.
  4. Evacuate the container according to the reaction needs. For the triple reaction vessel, the vacuum pump can be applied to evacuate the outer layer, to achieve the thermal insulation functionality.
  5. Set the temperature control device according to the temperature shown, to obtain the desired reaction temperature inside the container
  6. Slowly adjust the knob to achieve the desired rotation speed.
  7. Let the water or coolant pass through the condenser according to the process requirements.
  8. After the reaction, first turn the heating or cooling temperature control device function off, maintain the cycle and then close the circulation function when it is near to the ambient temperature.
- \*Note:** The temperature control unit cycle is turned off. The circulation valve cannot be closed, leave it open. Otherwise, if the body of the container is too hot or too cold, it could cause inflation and cause damages to the equipment.
9. After the speed test is set to 0, click on the key 1 minute after closing the reactor power
  10. Open the discharge valve to discharge the materials. If necessary, the valve parts or the sealing parts of the container should be cleaned and then reassembled safely.

**\*Note:** The glass parts must be cleaned after returning to normal temperature.

The droplets on the inside and outside should be cleaned after washing to avoid rupture of the container, caused by an irregular temperature or by a sudden change in temperature.

### 3. Technical Data

Table 3 - RE-650 speed-regulation glass reactor technical data

Model	RE-650
Stirring speed (rpm)	20~500
Jacket pressure (Mpa)	Inlet<=0.03
Pressure in the vessel (Mpa)	-0.1~0.0
Tolerable temperature difference (°C)	-80~200
Motor rated power (W)	370
Glass vessel capacity (L)	50
Jacket capacity (L)	About 3
Vessel jacket heat exchanging area (m <sup>2</sup> )	0.77
Condenser heat exchanging area (m <sup>2</sup> )	0.43
Diameter of circulating fluid inlet and outlet	DN15
Charging bottle (L)	2
Power supply	100 - 240V~, 50/60Hz
Ambient temperature (°C)	5~35
Relative humidity (%)	≤ 65
Protection degree	IP20
Dimensions (mm) WxDxH	900×690×2050
Drain port ground clearance (mm)	About 320
Net weight / Gross weight (kg)	82/124

#### 3.1 Borosilicate Glassware 3.3

- Chemical properties
- Hydrolytic resistance
- Acid resistance
- Alkali resistance
- Temperature resistance
- Optical

Technical features of working use in standard reactors: RE-650

Vacuum: -0.9 to +0.5 pressure bars

Working temperature: -30°C to +200°C

PTFE gaskets: -100°C a +250°C (maximum working -20 to +200°C)

Glass is an organic mixture of metal oxides fused together at high temperatures, which upon cooling, solidifies into the clear, rigid, non-crystalline, versatile material known widely across the globe

Borosilicate glass has a very high resistance to attack from water, acids, salt solutions, halogens and organic solvents. Only hydrofluoric acid, hot concentrated phosphoric acid and strong alkaline solutions cause appreciable corrosion of the glass.

Borosilicate glass has excellent thermal properties at both high and low temperatures. The maximum recommended working temperature for laboratory glassware manufactured is 500°C (for short periods of time only). Special care should be taken at temperatures above 150°C to ensure both heating and cooling is achieved in a slow and uniform manner.

## 3.2 Lubrication

High vacuum grease

Range Hg: 10-6 mm

Range: -40°C to 200°C invaluable

This grease does not melt and is been composed for lubrication and sealing of glass and wrenches in high vacuum systems with outstanding stability.

**Notice to customers:**

The product is made up of various components and various materials that must be recycled or, failing that, deposited in the corresponding debris removal sites when the product's life has been completed or when otherwise it is necessary to dispose of it. To do this, the end user who acquires the product must know the current regulations of each municipality and / or locality based on the waste electrical and electronic equipment. The user who acquires this product must be aware of and responsible for the potential effects of the components on the environment and human health as a result of the presence of hazardous substances. Never place the product in a conventional container of citizen scope if a previous dismantling and knowledge of the components that incorporates. If you do not know the procedure to follow, consult with the city council for more information.

## 4. Operating

### 4.1 Caution

Table 4.1 Caution

 Note	Please read this manual carefully before using it.
	Put it on the floor then lock the casters.
	The best ambient temperature for this device is 5~35°C. It will be broken if working outside this temperature range for a long period of time.
	Avoid strong vibration, otherwise it may damage the glass parts.
	Please connect the power supply according to the requirements on the name plate.
	Do not press the power cord. When there is a heavy thing on the power line or through the object, it may cause damage to the power cord and cause electric shock or fire.
	Please unplug the power plug from the socket to avoid long-term accumulation of dust if you do not use it for a long time, if the temperature is very high it may cause a fire.
	Do not pull the power cord strongly.
	Please disconnect the power supply and unplug the plug when moving it. Do not move it if there is power to avoid electric shock or fire.
	Please clean the glass parts before using it.
	When the glass device is connected, the interface should be uniformly applied to the vacuum grease to ensure the sealing effect.
	Please be careful when install the glass parts.
	Do not hit the glass.
	The viscosity of the circulating liquid must be less than 50cp when the reactor connected with dynamic temperature control system.
 Warning	Please start the circulating pump first, then start the Dynamic Temperature Control System.
	Do not heating the reactor with steam.
	Do not make the stirring paddle rotate at a high speed of no-load.
	Please contact us or searching for technical support when there's any abnormal situation.
 Warning	The maximum temperature difference between the jacket and the vessel is 80°C.
	Do not move the device when it is working.
	Be aware of scaling! Please wear glasses when dealing with the sample.
 Warning	Do not move it under high temperature!
	Do not refit this device.
	Do not use it outdoor. Get wet in the rain or splashing water on the surface may cause personnel injury.

## 4.2 Operating

Note:

Before starting the stirring function for the first time, pay attention to the relative position of the stirring paddle and the bottom of the vessel body to avoid collision.

### 4.2.1 Power Connection

When the reactor is connected to the power supply properly, the control panel will enter the welcome interface. 5 seconds later, it will change to the monitoring interface, which displays the current temperature, stirring speed and torque, as shown in the image.

### 4.2.2 Stirring speed and time setting

Press the Set knob on the control panel to enter the speed setting interface, turn the knob to the target value (clockwise to increase, counter clockwise to decrease) and then press again the Set knob to save the settings. Press it again to enter the time setting interface. Turn the knob slowly to set hour, minutes and seconds in turn, according to the actual needs. After each parameter is adjusted to the target value, they need to be saved by pressing the knob. If you don't press the Set knob, it will quit the current setting and return to the monitoring interface.

Note: When the reactor is empty, the stirring speed should be no more than 200rpm.

### 4.2.4 Operation

11. Feed liquid or solid materials into the reactor vessel slowly through the right port.
12. Click button Start/Stop on the control panel to start stirring speed control.
13. If the reaction should be carried on under vacuum conditions, start the vacuum pump to pump out the air in the vessel until it reaches vacuum state.
14. Start the Dynamic Temperature Control System according to actual needs of the reaction, and the temperature in the vessel can be controlled automatically when you set it ready.
15. When the reaction process is finished, shut down the Dynamic Temperature Control System when the temperature inside of the reactor is close to room temperature.

Note:

After shutting down the temperature control unit, please do not close the valves in the middle of the circulation hoses, keep all the hoses clear. Otherwise, when the temperature in the jacket or in the vessel changes dramatically, it may cause any crack or broken to the reactor body.

16. Shut down the vacuum pump.
17. Adjust stirring speed to minimum value; click button Start/Stop to stop stirring.
18. Click button Power On/Off to power off the controller.
19. Discharge the material from the draining port (14). After emptying the material, please disassemble the draining valve and clean the glass part and the PTFE stopper.
20. Reinstall the draining valve.

Note: When discharging the materials, slowly turn the black stopper at the bottom of the discharge valve counterclockwise. After cleaning the glass parts, wipe the liquid on the surface with a soft clean cloth to avoid contamination.

## **5. Maintenance and Management**

---

Ensure to make the equipment running normally to increase its service life, users should do daily maintenance and management. Before maintenance, please disconnect equipment power supply, preparing tools, materials according to according to the specific requirements in the operation manual. Otherwise, it will cause electric shock or damage.

### **5.1 Management**

Please follow following recommendations to keep the normal operation of the reactor:

- 1) Check the ambient temperature and humidity regularly to ensure a better working conditions.
- 2) Avoid severe vibration when moving it.
- 3) Keep the reactor from impact which might cause glass broken. Inspect the circulation hoses, replace the damaged or aged ones.

### **5.2 Cleaning**

- 1) Before cleaning the unit, disconnect the power cord. Otherwise, it may cause electric shock or fire hazard.
- 2) Use soft cloth to clean the surface.
- 3) Cleaning of glass parts should be done according to laboratory regulations.
- 4) Beware of hazardous substance leaks on the surface of the vessel or enters it.
- 5) Do not use brushes, grinding powder, acid, petrol or other solvents to clean the surface of the body to avoid damage to the protective layer.
- 6) The reactor body and components may be very hot, beware of burns. Clean it when the device glass is close to room temperature.
- 7) In order to avoid damage of components, please do not press the monitor and knob forcedly when cleaning them.

### **5.3 Maintenance**

The warranty of the equipment against manufacturing problems is of 12 months from the date of purchase. If the equipment has been used improperly, compensation for the damage caused will be charged.

Contact the Technical Service before obtaining professional maintenance to return the equipment to the factory:

1. Clean the machine so as not to damage the maintenance personnel.
2. Provide the description of the equipment failure.
3. Pack the machine properly to avoid damage during transport.

## 5.4 Troubleshooting

Table 5.3 Faults and troubleshooting

Faults	Reason	Troubleshooting
No display on touch screen after power on	Power supply is not connected	Check the power supply circuit
	* Fuse fault	Check the fuse and its installation
	Communication failure	Keep communication lines well connected
	PCB board fault	Please contact us or professional maintenance personnel
Displayed LAL on screen	Detected temperature is below the lower limit	Please contact us or professional maintenance personnel
Displayed HAL on screen	Detected temperature is above the upper limit	
Displayed OVER on screen	Temperature sensor failure	
Displayed ER-1 on screen	Module fault	
Displayed ER-2 on screen	Motor blocking fault	Inspect the voltage
Displayed ER-3 on screen	Hall logic error	
Displayed ER-4 on screen	Under voltage	
Displayed ER-5 on screen	Over voltage	
Displayed ER-6 on screen	Communication failure	Please contact us or professional maintenance personnel
Abnormal noise in stirring process	Stirring paddle installed improperly	Re-install stirring paddle
	Motor fault	Please contact us or professional maintenance personnel
	Debris in the vessel	Clean debris
Vacuum degree decrease	Glass joints are not well connected	Clean the joints and connect them properly
	Sealing ring wear	Replace sealing ring
	Sealing ring installed improperly	Reinstall sealing ring
	Vacuum hose aged	Replace vacuum hose
	Vacuum pumping capacity decreased	Replace vacuum pump
Discharge valve leak air	Dramatic temperature changes affected discharge valve sealing	Adjust the flange for sealing check
	Sealing ring damaged	Replace sealing ring

Note: \* The fuse is fixed above the power socket on the back of the motor controller. Two parts are combined as a whole unit.

## 6. Accessories and Spare Parts

Digital thermometer «TC-9226-A».  
Temperature range: from -200°C up to 1370°C  
**Part number 5903030**

Criohermostat "FRIGITERM-TFT-30"  
Accessory for the coolant circuit  
For regulable temperatures from -30 up to 100°C  
**Part number 3000547**

Gable lid  
Thermostatic bath accessory, which limits evaporation  
**Part number 3001295**

Thermostatic bath DIGITERM-TFT-200 -20  
Accessory for the reactor  
To regulate temperatures from ambient +5°C up to 200°C  
**Part number 3000550**

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Water recirculator  
"Interheater" Accessory of the  
coolant circuit  
To regulate temperature from +5° up to 90°C  
**Part number 6001491**

Rotary vacuum pump  
Flow volume 3.6 m<sup>3</sup>/h. Vacuum limit 0.1 mbar  
**Part number 5900620**  
spare oil for pump  
**Part number 5515030**

Vacuum membrane pump N-820  
Special for solvents  
Flow volume 20 litres/minute. Vacuum limit 8 mbar. PTFE rotors. PTFE coated  
membranes, valves in FPPM, motor protection level IP-44  
**Part number 5910009**

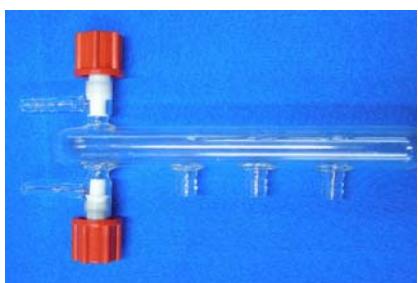
Scrubber unit  
Glass trap in Pyrex to avoid aspiration of liquids by means of the vacuum pump  
**Part number 4001611**

Universal clamp  
In dur-al with jaws in cork coated  
**Part number 7000403**

Double nut  
In chrome plated  
**Part number 7000408**

Double glass valve  
With three vacuum connection points (coolant, load & drip)  
**Part number 47078**

Glass T  
For vacuum connection  
**Part number 47077**

Silicone pipe  
External diameter 14 mm. Internal diameter 8 mm  
**Part number 46041**

Corrugated tube  
Reinforcement for the silicone tube  
**Part number 46079**

Fixing clamps  
Bags of 100 units  
**Part number 28006 (unit)**

PTFE bellows diameter 38\*40 x 75mm  
**Part number 1001984**

Conical connection cover 29/32 PTFE  
**Part number 28515**

Glass sensor adapter for reactor  
Combines with PH/T°C analyser part number 4120800  
& sensor part number 120801  
**Part number 4120802**

Guide for PTFE stirrers 29/32 Ø 10 MM  
High chemical resistance  
Low vacuum (5mmHg)  
Pressure (3-5 psi)  
High performance PTFE/PEEK/glass precision bearing seal  
Internal anti-whip sleeve  
**Part number 1001788**

Addition funnel 250ml PTFE key Step 4mm  
**Part number 1001795**

Load faucet 29/32 Step 4mm Olive 12/13  
**Part number 1001793**

Stainless steel funnel with handle 200 mm  
**Part number 1000909**

Support accessories trolley  
Total height 90 cm  
Lower platform dimensions: 70x70 cm  
Upper platform dimensions: 70x40 cm  
**Part number 1001744 (unit)**

Stainless steel bucket 15 litres  
**Part number 1001600**

Large opening Hoffman clamp  
**Part number 1000064**

50 litres reservoir  
**Part number 0106006**

#### OTHER ACCESSORIES:

- 0046030 PTFE tube 10X8 (1 meter)
- 0046149 Silicon tube to unload 15x20. (1 meter)
- 0046150 Silvar tube 14.4x8 mm reinforced with fiberglass (5 meters)
- 0046156 Silvar tube 19x12 mm reinforced with fiberglass (5 meters)
- 0046084 Viton tube 10x14 mm (1 meter)
- 0046085 Viton tube 12x17 mm (1 meter)
- 0046041 Silicon tube 14x8 mm (5 meters)
- 9515031 Fluid 47 V50 4 litres (silicone oil)
- 1001853 PTFE corrugated tube DN12 connection Ø10 (1 meter)
- 1001915 High vacuum grease 50 gr

## 7. Examples of Use of Support Trolleys

Frigiterm TFT & Interheater

Digiterm TFT & Interheater

Digiterm TFT, Vacuum pump, Scrubber & steel cubes  
Vacuum pump, Scrubber and 50 litres reservoir

## 8. Examples of Application